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## 30W Interchangeable AC Clip Adapter

PSAA30R-XXX-R Series

● 12V/15V/24V/56V ● DoE 6 ● Interchangeable AC Clip ● Cable ● Black

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### Features :

<b>Efficiency</b>	DoE Level VI (86.95% avg)
<b>Power Saving</b>	<0.1W @230Vac/50Hz
<b>Protection</b>	SCP, OVP, OCP
<b>Turn On Delay Time</b>	<3s @ 115Vac, max load
<b>Hold-up Time</b>	>10ms @ 115Vac, max load
<b>Rise Time</b>	<100ms @ 115Vac, max load
<b>Isolation</b>	Primary to Secondary 3000Vac,10mA for 1min.
<b>Insulation</b>	Primary to Secondary >7M ohm 500Vdc
<b>DC Output Cable (12V)</b>	1500±50mm, UL2468, AWG#18
<b>(15V, 24V, 56V)</b>	1500±50mm, UL2468, AWG#24
<b>DC Output Connector</b>	5.5x2.1x10mm, center positive, straight barrel
<b>Dimension</b>	82.7*55*39.1 (mm)
<b>Weight</b>	180 (g)

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declares that the following product  
 Product name: **PROFILER SCD**  
 Article no.: **00-76-900-4830**  
 Product type: **Digital readout unit**  
**for 2D and 3D measurements**

is in conformity with the following European Union harmonisation legislation, when operated in accordance with the operating manual:  
**EMC directive 2014/30/EU**  
**Low voltage directive 2014/35/EU**  
**RoHS directive 2011/65/EU**

The following harmonised standards have been applied:  
**DIN EN 61010-1:2020-03**  
**DIN EN 61326-1:2013-07**

In case of a modification of the product, which was not approved by the manufacturer, this Declaration of Conformity loses its validity.  
 This Declaration of Conformity is issued under the sole responsibility of the manufacturer.

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**Approved by** Uwe Schnitter | 26.04.2022  
**Issued by** Christopher Koch | 26.04.2022

**Proved to be QMS conform by** Uwe Schnitter | 26.04.2022  
**QM document** AB502\_QF07\_07 | Rev. A | Ver. 00 | Page 1 of 1

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Zubehör 10

## Table of Contents

1	Basics of the Operating Manual .....	109
1.1	Symbols and Abbreviations .....	110
2	Scope of Delivery .....	111
2.1	Optional Accessories .....	112
3	Product Description .....	113
3.1	Identification .....	113
3.2	Area of Application .....	113
3.3	Compliance with Directives and Standards .....	114
3.4	Structure and Ports .....	114
4	Safety .....	116
4.1	Intended Use .....	116
4.2	Installation Requirements .....	116
4.3	General Safety Instructions .....	117
4.4	Warning Note Categories .....	119
4.5	Hazard Symbols .....	119
5	Putting into Service .....	120
5.1	Putting the PROFILER SCD into Service .....	120
5.2	Connecting a 3rd Measuring System (Option) .....	122
5.3	Carrying out a Basic Configuration .....	125
5.4	Establishing the Connection to a PC .....	126
5.5	Connecting the Foot Switch (Option) .....	131
6	Operation .....	133
6.1	Becoming Familiar with the User Interface .....	133
6.2	Getting to Know the Position Display .....	134
6.3	Performing Measurements .....	137
6.4	Sending Position Values to a PC .....	138
6.5	Using the Position Correction .....	140
6.6	Enabling/Disabling the Standby Mode .....	151
6.7	Configuring the PROFILER SCD .....	152

7	Operation on a PC .....	171
7.1	GetPos Software .....	171
7.2	Instruction Set.....	180
8	Cleaning .....	181
9	Maintenance.....	182
10	Troubleshooting .....	183
11	Service .....	188
11.1	Replacement Components .....	189
12	Taking out of Service .....	190
13	Disposal .....	191
14	Technical Data.....	192
14.1	General Data.....	192
14.2	Encoder Interface.....	194
14.3	Ports.....	195
14.4	External 24 V/30 W Power Supply .....	202
15	Declaration of Conformity.....	204
16	Index.....	205

## 1 Basics of the Operating Manual

This operating manual gives you the information which is needed for trouble-free and safe use of the PROFILER SCD. You must have fully read and understood this operating manual and especially the safety instructions before you put into service, operate or dispose of the PROFILER SCD.

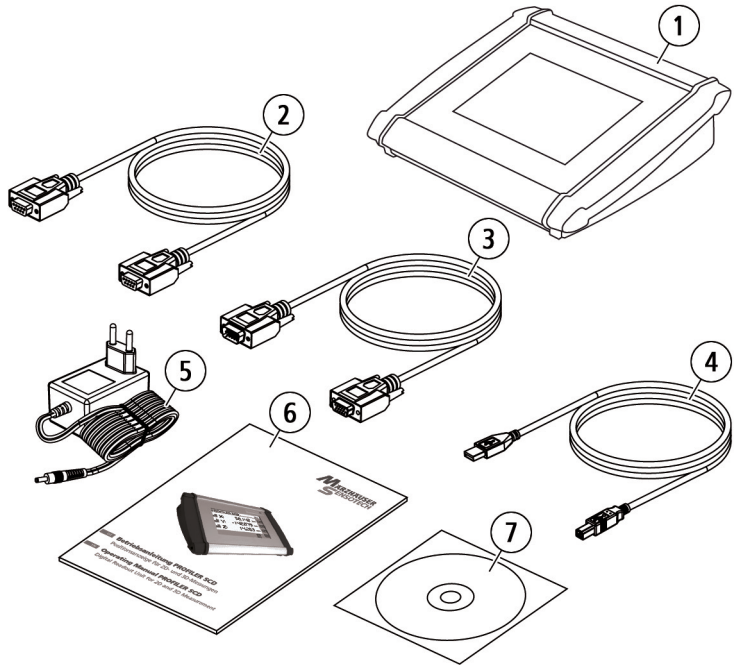
Märzhäuser Sensotech accepts no liability for damage and operating faults which result from ignoring this operating manual.

This operating manual is applicable for a PROFILER SCD with firmware version 1.26 and higher.

## 1.1 Symbols and Abbreviations

Symbol	Meaning
	additional information and useful notes
	start of an instruction
1, 2, 3 ...	steps within an instruction
»	cross reference to another part of this operating manual, another document or a website
Abbreviation	Meaning
approx.	approximately, roughly
e.g.	for example
fig.	figure
max.	maximum
min.	minimum
resp.	respectively
typ.	typical

## 2 Scope of Delivery

No.	Designation	Quantity
1	PROFILER SCD	1
2	RS-232 interface cable (9-pin)	1
3	MR encoder cable (15-pin, 2 channels, 2 m)	1
4	USB interface cable	1
5	External power supply 24 V/30 W (fig. similar) incl. 4 adapters (EU/GB/USA/AUS, no fig.)	1
6	Operating manual	1
7	Installation CD incl. drivers, operating manual, instruction set, software <i>GetPos</i>	1

© Mahaffey (Fig. 2), Samtec (Fig. 3), Kimblier/Walter (Fig. 4), source: 3D Content Central, Dassault Systèmes

## 2.1 Optional Accessories

The following optional accessories are available for the PROFILER SCD:

Designation	Order no.
Foot switch with 2 pedals	00-76-900-0862
Individual encoder cables for a 3rd measuring system	on request
Separate measuring systems (MR/analog 5Vss, 1Vss)	on request
Wall bracket	on request

For questions about ordering, please contact our sales team:

Tel.: **+49 6441 9116-10**

Fax: **+49 6441 9116-40**

E-Mail: **[sales@marzhauser.com](mailto:sales@marzhauser.com)**

If you want to make your own cable, you can find the pin assignments of all the PROFILER SCD ports here:

» *Chapter 14.3 Ports, page 195*

## 3 Product Description

### 3.1 Identification

You can find all required information for clear identification of the PROFILER SCD on the identification plate. The identification plate is attached to the underside of the PROFILER SCD.

### 3.2 Area of Application

The PROFILER SCD is used as a digital readout unit for measuring objects with a measuring stage from Märzhäuser Wetzlar.

Up to 3 incremental measuring systems can be connected to the PROFILER SCD. If needed, in addition to the X- and Y-axis of the measuring stage, an additional 3rd measuring system can be used.

Using the integrated position correction, systematic length errors of the measuring stage can be corrected. Evaluating reference signals creates an absolute reference of the position correction values to the machine zero.

The PROFILER SCD is operated via the integrated touchscreen. The PROFILER SCD can also be connected to a PC over an RS-232 or USB interface. The analysis and programming of the PROFILER SCD via the PC can be carried out with the *GetPos* software or directly by using the instruction set.

The PROFILER SCD is compatible with all measuring stages from Märzhäuser Wetzlar with an integrated measuring system.

Here you can find further information on the magnetic measuring system from Märzhäuser Sensotech:

» [www.marzhauser-st.com](http://www.marzhauser-st.com)

### 3.3 Compliance with Directives and Standards

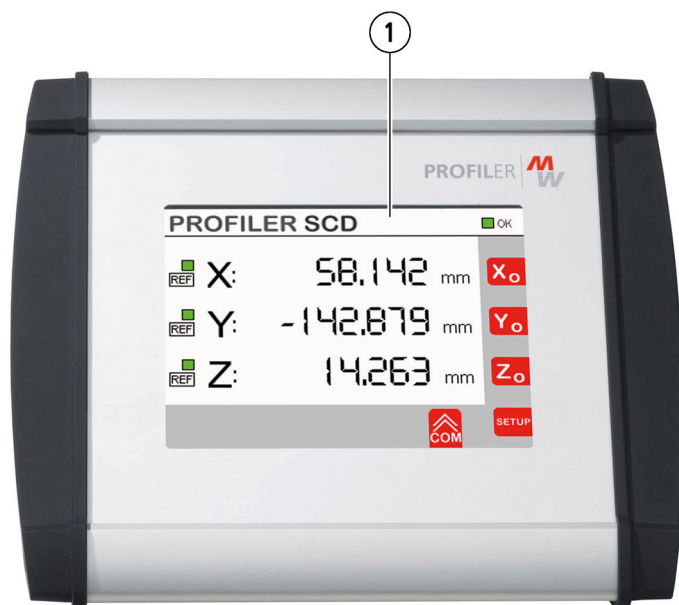
Märzhäuser Sensotech hereby declares that the PROFILER SCD was designed and produced in compliance with the EU directives and harmonized standards mentioned in the Declaration of Conformity.


The Declaration of Conformity can be found here:

» *Chapter 15 Declaration of Conformity, page 204*

### 3.4 Structure and Ports

No.	Designation	Function
1	-	touchscreen

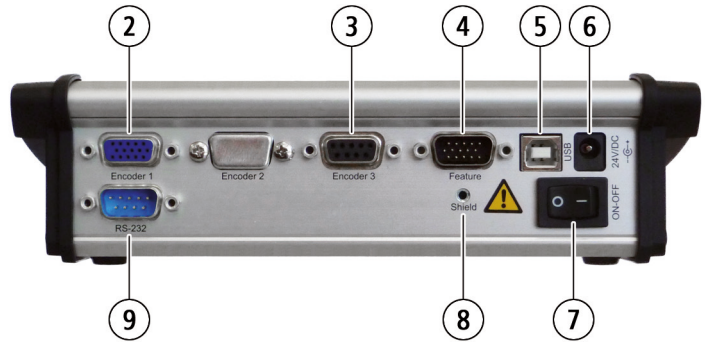No.	Designation	Function
2	Encoder 1	port for measuring stage with 2 measuring systems ( <i>MR/analog 5Vss</i> )
3	Encoder 3	port for 3rd measuring system ( <i>MR/analog 5Vss, 1Vss, TTL (RS-422)</i> )
4	Feature	optional functions (port for foot switch, reference switch input for 3rd measuring system)
5	USB	USB port (PC interface)
6	24V/DC	port for power supply (24 V)
7	ON-OFF	on/off switch
8	Shield	port for shield/ground
9	RS-232	RS-232 port (PC interface)

## 4 Safety

### 4.1 Intended Use

The PROFILER SCD is intended and designed for the connection and evaluation of

- 1 measuring stage from Märzhäuser Wetzlar with 2 incremental measuring systems type *MR/analog 5Vss* and
- 1 further incremental measuring system without external power supply type *MR/analog 5Vss, 1Vss* or *TTL (RS-422)*.

On request, your PROFILER SCD can also be upgraded for the connection of an incremental measuring system of type *TTL (non-differential)* on the port *Encoder 3*. When needed, please contact our Customer Service Center:

» *Chapter 11 Service, page 188*

### 4.2 Installation Requirements

Category	Condition
Installation location	closed rooms, no risk of explosion
Installation height	max. 2,000 m
Environmental temperature	+5...+40 °C
Storage temperature	-20...+70 °C
Cooling	normal convection
Humidity	max. 85%, non-condensing

## 4.3 General Safety Instructions

- Only use the PROFILER SCD for the purposes which are described in this operating manual.
- Read and understand the operating manuals of the measuring systems and peripherals, which you would like to connect to the PROFILER SCD.
- Only connect measuring stages and measuring systems from Märzhäuser Wetzlar to the port *Encoder 1*.
- Only connect incremental measuring systems of type *MRI analog 5Vss, 1Vss or TTL (RS-422)* without external power supply to the port *Encoder 3*.
- Only use shielded cables for connecting measuring systems, PCs and other peripherals.
- The PROFILER SCD is insulated to the power supply. Ensure that no dangerous voltages or currents are induced by peripheral devices in the PROFILER SCD.
- Only use the supplied 24 V/30 W power supply to power the PROFILER SCD.
- Ensure that the mains voltage matches the label on the power supply before connecting the power supply to the mains.
- Disconnect the PROFILER SCD from the mains before connecting/disconnecting measuring systems, a PC or other peripherals to/from the PROFILER SCD.
- Set the PROFILER SCD in such a way that you can immediately disconnect the power supply from the mains, without hindrance, in case of emergency.
- Ensure that no chips, liquids, objects or substances enter the interior of the PROFILER SCD or the power supply. Protect the PROFILER SCD from dust and dirt as much as possible.

- The PROFILER SCD is not protected against corrosive, infectious, radioactive or other substances, which are harmful to health. When handling such substances, ensure that all legal requirements are fulfilled, especially the national regulations on accident prevention.
- Immediately replace the power supply if there is damage on the housing or cable of the power supply.
- Disconnect the PROFILER SCD from the mains if there are malfunctions, defects or safety flaws. Secure the PROFILER SCD against restarting.
- Never open the housing of the PROFILER SCD or the accompanying power supply. In case of violation, all warranty and liability claims will become void.
- Only have repair work carried out by our Customer Service Center.  
» *Chapter 11 Service, page 188*
- Ensure that the PROFILER SCD, in conjunction with your application, meets the applicable safety regulations and legal requirements.
- Keep this operating manual to hand in the place in which the PROFILER SCD is used.
- If you lend or sell the PROFILER SCD, you must also provide all the documents supplied with it.

## 4.4 Warning Note Categories

Warning notes are divided into the following categories:

### **WARNING**

Hazardous situation that can cause serious injury or death if not avoided

---

### **NOTICE**

Warning of possible damage or malfunction

---

## 4.5 Hazard Symbols

The following hazard symbols are used on the PROFILER SCD or in warning notes in this operating manual:

<b>Symbol</b>	<b>Meaning</b>
	warning of a general hazard, which can be avoided by observing the operating manual
	warning of hazard from electric current
	warning of hazard from substances which are harmful to health

## 5 Putting into Service

### 5.1 Putting the PROFILER SCD into Service

#### **NOTICE**

**Possible equipment damage from putting into service / taking out of service with existing power connection.**

Disconnect the PROFILER SCD from the mains before connecting/disconnecting measuring systems, a PC or other peripherals to/from the PROFILER SCD.

**Possible incorrect data transmission from using unshielded cables.**

Only use shielded cables for connecting measuring systems, PCs and other peripherals.

---

#### ► Putting the PROFILER SCD into service

- 1 Read and understand the *Safety* chapter before installation and putting into service.  
» *Chapter 4 Safety, page 116*
- 2 Read and understand the operating manuals of the peripherals, which you would like to connect to the PROFILER SCD.
- 3 Assemble your measuring stage in accordance with the Assembly Instructions.
- 4 Ensure that the PROFILER SCD has a secure footing and is assembled stably.
- 5 Ensure that the PROFILER SCD is switched off (switch *ON-OFF* to position *0*).

**NOTICE****Possible equipment damage from connecting inappropriate measuring systems.**

Only connect measuring stages and measuring systems from Märzhäuser Wetzlar to the port *Encoder 1*.

- 6 Connect your measuring stage to the port *Encoder 1* of the PROFILER SCD. Use the MR encoder cable for this.

**Note:** The reference switches of your measuring stage are thus also connected to the PROFILER SCD.

**WARNING****Possible electric shock and equipment damage from connecting foreign or damaged power supplies.**

Only use the delivered 24 V/30 W power supply. Immediately replace the power supply if there is damage on the housing or cable of the power supply. Only use replacement components from Märzhäuser Sensotech.

**Possible electric shock and equipment damage from connecting to non-specified mains.**

Ensure that your mains voltage matches the label on the power supply before connecting the power supply to the mains.

- 7 Connect the power supply cable to the port *24VDC*.
- 8 Connect the power supply to the mains.
- 9 Switch on the PROFILER SCD (switch *ON-OFF* to position *I*).

The display lights up. The PROFILER SCD starts the boot sequence and is operational after 1–2 seconds.

- 10** If you have received the PROFILER SCD as an accessory for a measuring stage, the PROFILER SCD is already specifically pre-configured for your measuring stage.  
If you have received the PROFILER SCD independently from a measuring stage, carry out the basic configuration.
- » *Chapter 5.3 Carrying out a Basic Configuration, page 125*
- 11** Move the measuring stage at least 1 cm in both directions.
- Note:** The measuring systems are automatically calibrated in this way. The calibration is kept after switching off the PROFILER SCD and is always fed into the operation.

Putting into service is completed. You can now perform measurements with the PROFILER SCD.

If needed, adapt the PROFILER SCD to your individual needs:  
» *Chapter 6.7 Configuring the PROFILER SCD, page 152*

## 5.2 Connecting a 3rd Measuring System (Option)

In addition to your measuring stage, with the PROFILER SCD you can evaluate a 3rd measuring system (probe, linear encoder or rotary encoder).

On the port *Encoder 3*, you can connect an incremental measuring system without external power supply from Märzhäuser Sensotech (*MR/analog 5Vss*) or from another manufacturer with an interface of type *1Vss* and *TTL (RS-422)*.

You can receive the appropriate encoder cable for your 3rd measuring system as an optional accessory at our Customer Service Center:  
» *Chapter 11 Service, page 188*

### ► Connecting a 3rd measuring system

- 1 Read and understand the operating manual of your 3rd measuring system.

#### **NOTICE**

##### **Possible equipment damage from putting into service / taking out of service with existing power connection.**

Disconnect the PROFILER SCD from the mains before connecting/disconnecting measuring systems, a PC or other peripherals to/from the PROFILER SCD.

- 2 Switch off the PROFILER SCD (switch *ON-OFF* to position *0*).
- 3 Disconnect the power supply of the PROFILER SCD from the mains.

#### **NOTICE**

##### **Possible equipment damage from connecting inappropriate measuring systems.**

Only connect incremental measuring systems of type *MR/ analog 5Vss, 1Vss* or *TTL (RS-422)* without external power supply to the port *Encoder 3*.

##### **Possible incorrect data transmission from using unshielded cables.**

Only use shielded cables for connecting measuring systems, PCs and other peripherals.

- 4 Connect your 3rd measuring system to the port *Encoder 3* of the PROFILER SCD. Use an appropriate encoder cable for this.
- 5 If you use a reference switch in combination with your 3rd measuring system, connect the reference switch to the port *Feature*. Use an appropriate cable for this.
  - » *Chapter 2.1 Optional Accessories, page 112*

**Note:** If your 3rd measuring system has reference marks, the reference signal is transmitted over the port *Encoder 3* to the PROFILER SCD.

- 6 Connect the power supply of the PROFILER SCD to the mains.
- 7 Switch on the PROFILER SCD (switch *ON-OFF* to position *I*).

The display lights up. The PROFILER SCD starts the boot sequence and is operational after 1–2 seconds.

- 8 For the 3rd measuring system, carry out a basic configuration.
  - » *Chapter 5.3 Carrying out a Basic Configuration, page 125*
- 9 If you are using a measuring system of type *MR/analog 5Vss*, you must calibrate your 3rd measuring system. For this purpose, move your 3rd axis at least 1 cm in both directions.

**Note:** The calibration is kept after switching off the PROFILER SCD and is always fed into the operation. Measuring systems of type *1Vss* and *TTL (RS-422)* do not have to be calibrated.

Your 3rd measuring system can now be assessed by the PROFILER SCD.

## 5.3 Carrying out a Basic Configuration

If you have received the PROFILER SCD as an accessory for a measuring stage, and are not using a 3rd measuring system, no basic configuration is needed.

If, however, you have upgraded the PROFILER SCD or are using a 3rd measuring system, you must carry out a basic configuration for all measuring systems, which are not part of a measuring stage within the scope of delivery.

### ► Carrying out a basic configuration

- 1 Open the menu *Setup*.
- 2 In the menu *Encoder*, configure all measuring systems which are not part of a measuring stage within the scope of delivery.  
» *Chapter 6.7.6 Encoder, page 163*
- 3 In the menu *Origin*, configure all measuring systems which are not part of a measuring stage within the scope of delivery.  
» *Chapter 6.7.3 Origin, page 154*
- 4 If needed, perform a position correction for all measuring systems, which are not part of a measuring stage within the scope of delivery.  
» *Chapter 6.7.5 PosCorr (Position Correction), page 161*

The basic configuration of the PROFILER SCD is now completed.

## 5.4 Establishing the Connection to a PC

As an option, you can connect the PROFILER SCD to a PC over the RS-232 or USB interface. The interface is automatically switched in the PROFILER SCD.

If the PROFILER SCD is connected to a PC over the USB interface, the RS-232 interface is automatically disabled. Simultaneous use of both interfaces is not possible.

### 5.4.1 Establishing a USB Connection

#### WARNING

**Possible electric shock and equipment damage from connecting inappropriate peripherals.**

The PROFILER SCD is insulated to the power supply. Ensure that no dangerous voltages or currents are induced by peripheral devices in the PROFILER SCD.

---

With a USB connection, communication between the PROFILER SCD and the PC takes place via a virtual COM interface. To connect the PROFILER SCD with a USB cable to your PC, you need *Windows 7 (32/64 bit)*, *Windows 8 (32/64 bit)*, *Windows 10 (32/64 bit)* or *Windows 11 (64 bit)*.

Observe the following notes, regardless of the version of your operating system:

- Allow the drivers to install automatically from the operating system.
- Do not connect to *Windows Update*.
- Continue installing the drivers despite the message about the missing *Windows-Logo-Test*.

► **Establishing the connection to a PC using a USB cable**

- 1 Switch off the PROFILER SCD (switch *ON-OFF* to position *0*).

## NOTICE

**Possible equipment damage from putting into service / taking out of service with existing power connection.**

Disconnect the PROFILER SCD from the mains before connecting/disconnecting measuring systems, a PC or other peripherals to/from the PROFILER SCD.

- 2 Disconnect the power supply of the PROFILER SCD from the mains.
- 3 Switch on the PC.
- 4 Insert the installation CD in the CD drive.
- 5 Connect the port *USB* of the PROFILER SCD with a USB port of your PC. Use the USB interface cable for this.
- 6 Connect the power supply of the PROFILER SCD to the mains.
- 7 Switch on the PROFILER SCD (switch *ON-OFF* to position *I*).

The operating system automatically recognises a new device (*PROFILER SCD*).

- 8 Follow the installation instructions of your operating system.

**Note:** If the driver for the PROFILER SCD is not automatically found on the CD, enter the archive location of the driver (.inf file). You can find the drivers for the PROFILER SCD on the installation CD in the *..Drivers* folder.

Once installation completes, the PROFILER SCD is connected to your PC.

Here you can read how you can use the PROFILER SCD together with your PC:

» *Chapter 7 Operation on a PC, page 171*

### ► Checking the version of the installed driver

If needed, you can check the version of the installed driver.

- 1 Open the Device Manager:
  - **Windows 7:** *Start → Control Panel → System → Device Manager*
  - **Windows 8:** right-click in the left, lower corner of the desktop/start screen → *Device Manager*
  - **Windows 10/11:** right-click on the Windows symbol in the left, lower corner of the desktop/start screen → *Device Manager*
- 2 Expand the column *Ports (COM and LPT)* by double-clicking.
- 3 Double-click on the entry *PROFILER SCD* to open the context menu.
- 4 In the context menu, select the entry *Properties*.
- 5 In the *Properties* window, select the tab *Drivers*.

In the tab *Drivers*, you can see the version of the installed driver.

## 5.4.2 Establishing an RS-232 Connection

### WARNING

#### **Possible electric shock and equipment damage from connecting inappropriate peripherals.**

The PROFILER SCD is insulated to the power supply. Ensure that no dangerous voltages or currents are induced by peripheral devices in the PROFILER SCD.

#### ► **Establishing the connection to a PC using an RS-232 cable**

- 1 Switch off the PROFILER SCD (switch *ON-OFF* to position *0*).

### NOTICE

#### **Possible equipment damage from putting into service / taking out of service with existing power connection.**

Disconnect the PROFILER SCD from the mains before connecting/disconnecting measuring systems, a PC or other peripherals to/from the PROFILER SCD.

- 2 Disconnect the power supply of the PROFILER SCD from the mains.
- 3 Connect the port *RS-232* of the PROFILER SCD with a serial port of your PC. Use the RS-232 interface cable for this.
- 4 Switch on the PC.
- 5 Connect the power supply of the PROFILER SCD to the mains.
- 6 Switch on the PROFILER SCD (switch *ON-OFF* to position *1*).

The PROFILER SCD is now connected with your PC.  
It is not necessary to install drivers with an RS-232 connection.

If you would like to read the PROFILER SCD over a terminal or other software, use the following settings to configure the COM interface:

<b>Parameter</b>	<b>Value</b>
Baud rate	57,600 baud
Data bits	8
Stop bits	1
Parity	none
Flow control	none

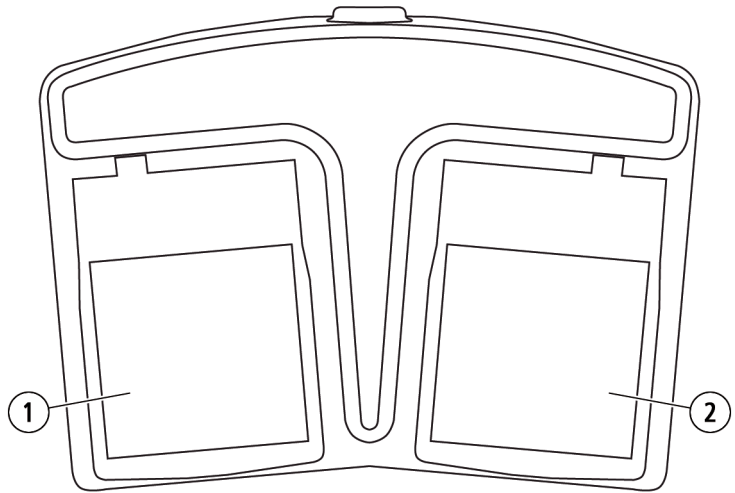
Here you can read how you can use the PROFILER SCD together with your PC:

» *Chapter 7 Operation on a PC, page 171*

## 5.5 Connecting the Foot Switch (Option)

The foot switch from Märzhäuser Sensotech enables hands-free operation of the PROFILER SCD. Using the foot switch, you can send the position values to a connected PC and zero position values.

No.	Designation	Function
1	left pedal	zero position values of all axes
2	right pedal	send position values to the PC

The foot switch is available as an optional accessory.  
 » Chapter 2.1 Optional Accessories, page 112

► **Connecting the foot switch**

- 1 Switch off the PROFILER SCD (switch *ON-OFF* to position *0*).

**NOTICE**

**Possible equipment damage from putting into service / taking out of service with existing power connection.**

Disconnect the PROFILER SCD from the mains before connecting/disconnecting measuring systems, a PC or other peripherals to/from the PROFILER SCD.

- 
- 2 Disconnect the power supply of the PROFILER SCD from the mains.
  - 3 Connect the foot switch cable to the port *Feature* on the PROFILER SCD.
  - 4 Connect the power supply of the PROFILER SCD to the mains.
  - 5 Switch on the PROFILER SCD (switch *ON-OFF* to position *1*).

The foot switch is now ready for operation.

## 6 Operation

### 6.1 Becoming Familiar with the User Interface

In manual mode, operate the PROFILER SCD over the user interface on the touchscreen.

Alternative operating options:


» Chapter 7 Operation on a PC, page 171

selection box    inactive check box    activated check box    deactivated check box

The screenshot shows the 'Encoder 1/2' configuration screen. It is divided into three columns for X, Y, and Z axes. Each column has a title box (X, Y, Z), an ON/OFF checkbox, a Count +/- checkbox, an Encoder type section with MR, TTL, and 1Vss checkboxes, a REF and Corr checkbox, and a Period [mm] input field. The X and Y axes have a Period of 0.500000, while the Z axis has 0.00025. Below the input fields is a numeric keypad with buttons 0-9, a decimal point, and a carriage return key. To the right of the keypad is a vertical bar with four buttons: a red SAVE button, a grey up arrow button, a red down arrow button, and a red ESC button. Labels on the right side identify these as 'active button', 'inactive button', and 'entry bar'. Labels at the bottom identify the input fields as 'inactive input field' and 'active input field'.


The *entry bar* appears as soon as you tap on an *active input field*. The following buttons are displayed in the *entry bar*:

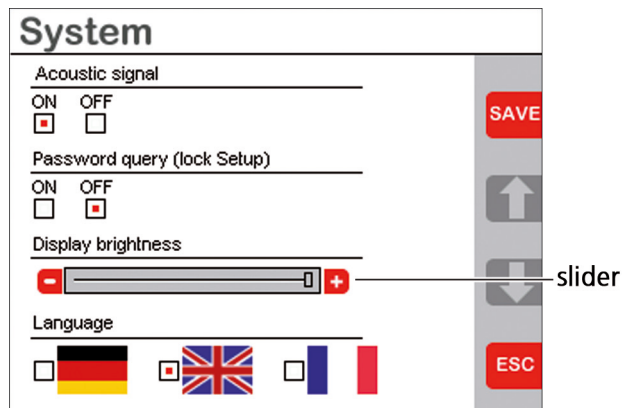-  enter numbers
  -  enter point
  -  delete last character (number or point)
  -  change signs
- Note:** The button  changes to  when all characters in an input field have been deleted.
-  confirm entry

## 6.2 Getting to Know the Position Display

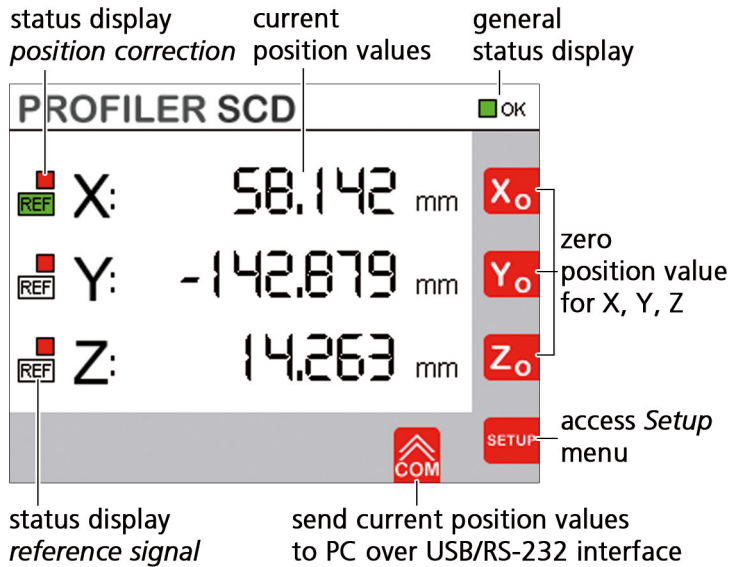The position display appears after switching on the PROFILER SCD. Which displays and buttons are shown in the position display depends on how your PROFILER SCD is configured.

The X- and the Y-axis of the measuring stage, which is connected to the port *Encoder 1*, are assigned to the position displays X and Y. If you have connected a 3rd measuring system to the port *Encoder 3*, the 3rd measuring system is assigned to the position display Z.

## 6.2.1 Current Position Values

The position display shows you the current position values of the measuring system with respect to the most recent zero point.

At the factory, the number of decimal places is pre-configured to 3 and the measuring unit to *mm*. You can, however, individually adapt the number of decimal places and the measuring unit to your needs.

» Chapter 6.7.4 Display, page 160

## 6.2.2 General Status Display

The general status display in the right-hand, upper corner of the touchscreen gives you information about the PROFILER SCD's readiness for operation.

 **OK Green**      The PROFILER SCD is ready for operation.

 **E4 Red**      The PROFILER SCD is not ready for operation. An error code is displayed next to the red display field.

Should the general status display be red, contact our Customer Service Center and give the displayed error code.

» *Chapter 11 Service, page 188*

## 6.2.3 Status Display Position Correction

If position correction was enabled in the menu *Encoder*, the display field  appears to the left of the respective position display (X, Y, Z).

 **Red, flashing**      Machine zero was not yet set or approached.  
The position correction is not active.

 **Green, constant**      Machine zero was set/approached.  
The position correction is active.

You can select the function for setting the machine zero in the menu *Origin*.

» *Chapter 6.7.3 Origin, page 154*

## 6.2.4 Status Display Reference Signal

If reference signal evaluation was enabled in the menu *Encoder*, the display field **REF** appears to the left of the respective position display (X, Y, Z).

<b>REF</b> White	The reference signal evaluation is enabled.
<b>REF</b> Green	A reference switch or reference mark was approached.

## 6.3 Performing Measurements

### NOTICE

#### Possible malfunction from incomplete putting into service.

Before you can take measurements with the PROFILER SCD, you must have fully put the PROFILER SCD into service.

» *Chapter 5 Putting into Service, page 120*

#### ► Performing measurements

- 1 If the position display shows *REF SEARCH*, proceed according to the function set for setting the machine zero (*Chapter 6.7.3 Origin, page 154*):
  - Move to the reference switches/reference marks of all axes or
  - set the machine zero with **X<sub>0</sub>**, **Y<sub>0</sub>** and **Z<sub>0</sub>**.
- 2 If the position correction is configured for your measuring stage, enable the position correction.
  - » *Chapter 6.5 Using the Position Correction, page 140*

- 3 Position the sample/workpiece below the measuring equipment (e.g. microscope) by moving your measuring stage.
- 4 Zero the position values with ,  or .
 

**Note:** With a connected foot switch (optional accessory), you can also zero the position values by pressing the left pedal.
- 5 Approach the desired measuring point.

Now you can read the current position values from the position display.

## 6.4 Sending Position Values to a PC

You can send position values to a connected PC in two different ways.

### ► Sending position values to a PC by a trigger event

- 1 Send the current position values to a connected PC via
  - button  on PROFILER SCD
    - » *Chapter 6.7.7 COM (Data / Interface), page 167*
  - foot switch (right pedal)
    - » *Chapter 5.5 Connecting the Foot Switch (Option), page 131*

The current position values are sent to a connected PC and can be received by GetPos, a terminal or an individual software solution.

### ► Sending position values to a PC by a position query

- 1 Send the current position values to a connected PC via
  - button  in GetPos
    - » *Chapter 7.1 GetPos Software, page 171*
  - button *Enter* on your PC keyboard while GetPos is running
    - » *Chapter 7.1.4 Key Layout on a PC, page 178*

- instruction
  - » *Chapter 7.2 Instruction Set, page 180*

The current position values are sent to a connected PC and can be received by GetPos, a terminal or an individual software solution.

On the PC, you can differentiate the trigger event from a position query by evaluating the first character of the sent information:

- **Example of a return value for a position query initiated by a trigger event:**

*>X 123.456 mm CR LF*

First character: > (0x3E hex)

Answer length: 16 characters<sup>1</sup> + CR LF

- **Example of a return value for a position query initiated on the PC:**

- **Instruction Set SCD<sup>plus</sup>**

*X 123.456 mm CR LF*

First character: Space (0x20 hex)

Answer length: 16 characters<sup>1</sup> + CR LF

- **Instruction Set PROFILER**

*2.500*

First character: Zahl

Answer length: variable

<sup>1</sup> If you have selected *mil* on the PROFILER SCD as the measuring unit to display the position value, a return value consists of 17 characters + CR LF.

## 6.5 Using the Position Correction

### **NOTICE**

#### **Loss of preconfigured position correction values from deleting or overwriting.**

The menu *Position Correction* is not write-protected. Do not delete, overwrite or expand any values entered by Märzhäuser Wetzlar. If required, protect the menu *Setup* using a password against unauthorised access.

» *Chapter 6.7.2 System, page 153*

#### **Incorrect position values from relocating the machine zero or changing the encoder type.**

Position correction values are valid as long as the location of the machine zero and the chosen encoder type are unchanged. If you, after entering position correction values,

- relocate the machine zero,
- mechanically offset a reference switch,
- move to another reference mark to set the machine zero,
- change the function for determining the machine zero or
- change the encoder type,

you must delete all previously stored position correction values and enter new position correction values.

---

### 6.5.1 Functioning of the Position Correction

The PROFILER SCD is equipped with a piecewise linear position correction, in order to compensate for length measurement errors of your measuring stage and an additional axis. The measurements can be corrected using up to 128 reference positions for each axis. The PROFILER SCD calculates position values between the reference positions through linear interpolation.

### 6.5.2 Using the Pre-Configured Position Correction

If you have received the PROFILER SCD as an accessory for a manual measuring stage with an integrated measuring system, the position correction of the PROFILER SCD is already configured for the X- and Y-axis of your measuring stage.

You can carry out measurements as soon as the machine zero is approached/set for all axes.

» *Chapter 6.5.6 Enabling the Position Correction, page 150*

### 6.5.3 Configuring the Position Correction Yourself

You need to configure the position correction by yourself if you

- have received the PROFILER SCD as an accessory for a measuring stage from Märzhäuser Wetzlar, but would like to connect an additional measuring system, or
- have purchased the PROFILER SCD to upgrade a measuring stage from Märzhäuser Wetzlar or
- want to use the PROFILER SCD with a measuring setup which is not manufactured by Märzhäuser Wetzlar.

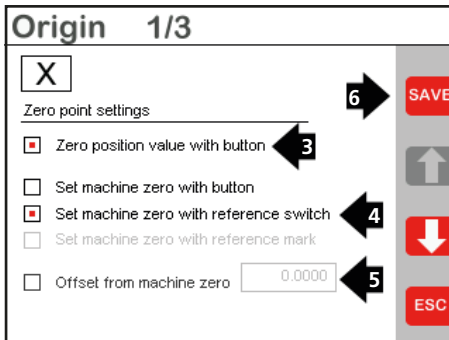
### Quick Overview

- A Configure the *Origin* settings.
- B Configure the *Encoder* settings. At first, it is necessary to disable the position correction.
- C Configure the position correction for all axes which need to be corrected one after another. For this, use a measurement scale placed on the measuring stage/setup. Feed the PROFILER SCD with value pairs of target positions and actual positions at discrete measuring points while moving the measuring stage/setup along the axes which need to be corrected.
- D Enable the position correction.
- E Restart the PROFILER SCD.
- F Set or move to the machine zero of your measuring stage/setup.
- G Carry out corrected measurements.


Now follow all of the following steps carefully.

► **Configuring the position correction yourself**



As an example, the position correction is performed for the X-axis within the following steps.

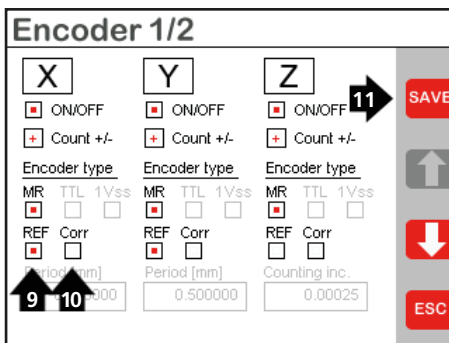- 1 Open the menu *Setup*.
- 2 Open the menu *Origin 1/3*.
- 3 Enable the function *Zero position value with button*.
- 4 Enable the desired function for setting the machine zero (Märzhäuser measuring stage: *Set machine zero with reference switch*).
- 5 Make sure that the function *Offset from machine zero* is disabled.

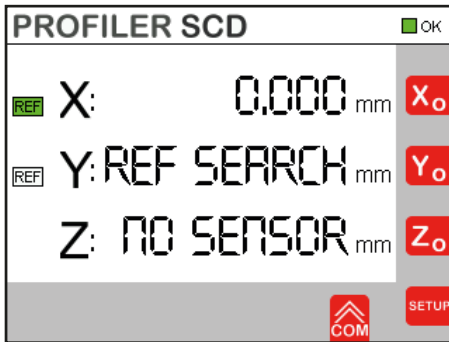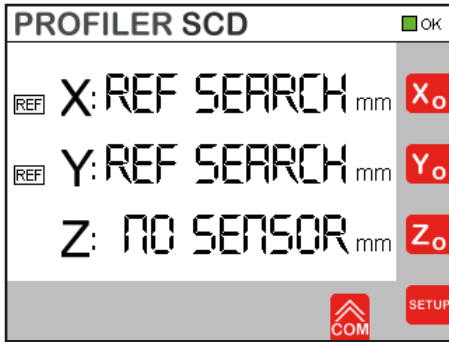- 6 Save your entries with **SAVE**.
- 7 Quit the menu *Origin 1/3* with **ESC**.
- 8 Open the menu *Encoder 1/2*.
- 9 Enable the *REF* check box.
- 10 Disable the *Corr* check box.
- 11 Save your entries with **SAVE**.
- 12 Quit the menu *Encoder 1/2* with **ESC**.
- 13 Quit the menu *Setup* with **ESC** to switch to the position display.

14 Place a measurement scale (e.g. glass scale) on your measuring stage/setup along the X-axis.

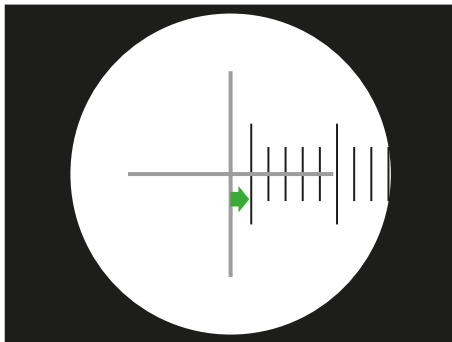
15 Move to the machine zero of the X-axis.

- If you have enabled in the *Origin* menu the function *Set machine zero with reference switch*, or *Set machine zero with reference mark*, move to the the reference switch or the desired reference mark of the X-axis.
- If you have enabled in the *Origin* menu the function *Set machine zero with button*, position your measuring stage/setup as desired. Set the machine zero with **X<sub>0</sub>**.


If the *status display reference signal* **REF** to the left of the X-axis' position display changes from white to green, the machine zero was successfully approached.

The position display of the X-axis switches from *REF SEARCH* to *0* (unit of measurements and number of decimal places depend on your individual settings).

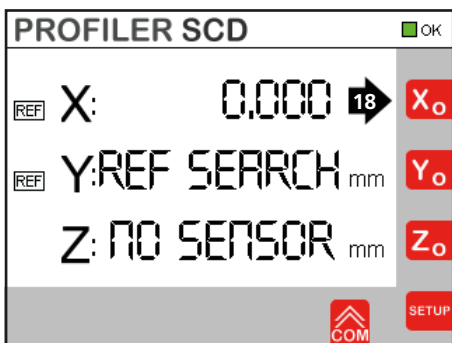
- 16 Make sure that the measurement scale is aligned parallel to the X-axis and that the scale of your measurement scale is visible through the optics of your microscope.
- 17 Move to the nearest measuring point available on your measurement scale.

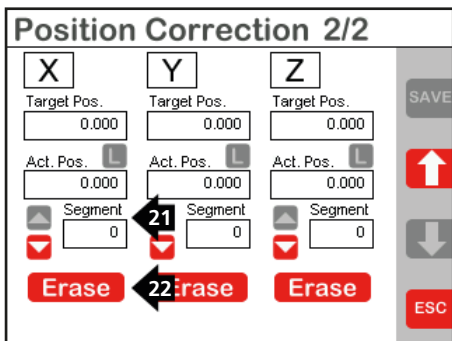
- 18 Zero the position value of the X-axis with **Xo**.

- 19 Open the menu *Setup*.
- 20 Open the menu *Position Correction (PosCorr)* and switch to the menu *Position Correction 2/2* with **↓**.
- 21 Delete all segments along with the associated target and actual values which are saved for the X-axis. To do so, navigate with **↑** to segment 0.
- 22 Delete all segments which are saved for the X-axis with **Erase**.
- 23 Confirm the query *Delete all position correction data of axis?* with **Yes**.

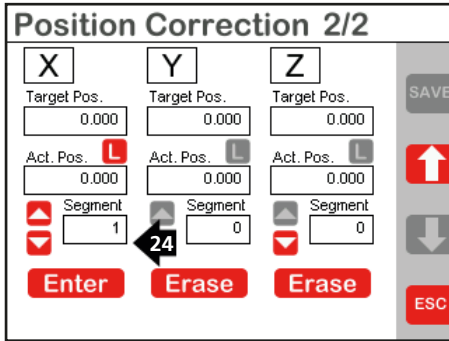
All segments which were saved for the X-axis, along with their associated target and actual values, are now deleted.

24 Add a new segment with .


**Note:** Target and actual value for the 0 segment are pre-defined and cannot be changed.

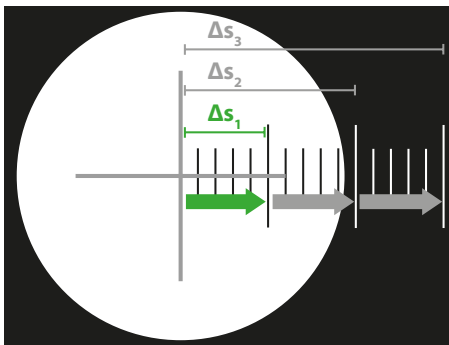The number of the new segment appears in the *Segment* display field.

25 Move the measuring stage/setup to the next measuring point of the measurement scale.

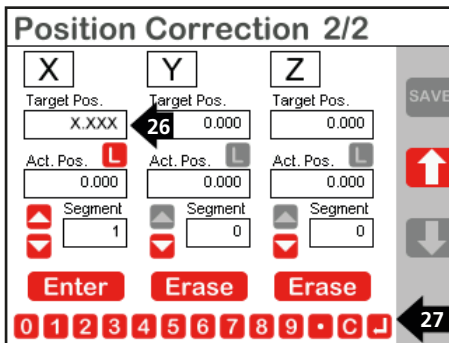**Note:** Select a constant distance between the measuring points depending on the travel range of your measuring stage/setup and your measurement scale. Example: With a travel range of 130 mm, a suitable constant distance between the measuring points can be 10 mm or 1/2 inch.

26 Tap on the *Target Pos.* input field.

27 Enter the distance  $\Delta s$ , travelled up to the current measuring point on the measurement scale, via the entry bar (for 1<sup>st</sup> segment:  $\Delta s_1$ , for all following segments:  $\Delta s_2$ ,  $\Delta s_3$ ,  $\Delta s_4$  and so on up to  $\Delta s_n$ ).

28 Confirm your entry with .


**NOTICE**

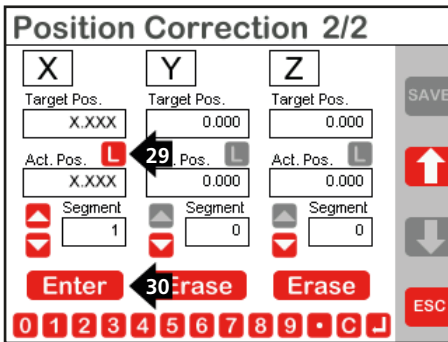**Incorrect position correction from excessive deviation between target value and actual value.**


The *Target Pos.* and *Act. Pos.* values may deviate from each other by a maximum of 127 µm. Check each value pair before saving it.

**Incorrect position correction from entering negative values.**

Only positive values may be entered for *Target Pos.* and *Act. Pos.* Do not enter values with a negative sign.

- 29 Adopt the currently measured position value in the input field *Act. Pos.* with **L**.
- 30 Save the entered target and actual value for the segment with **Enter**.
- 31 To define further segments (max. 128), repeat steps 24 to 30.
- 32 Quit the menu *Position Correction 2/2* with **ESC**.
- 33 To configure the position correction of the Y-axis and the Z-axis (if applicable), carry out steps 1 to 32 for the Y-axis and the Z-axis.
- 34 Remove the measurement scale from the measuring stage/setup.

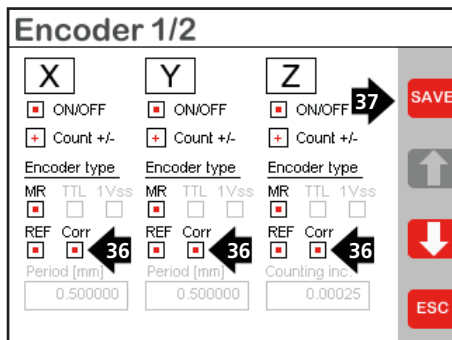
35 Open the menu *Encoder 1/2*.

36 Enable the check box *Corr* for the axes you have configured the position correction for.

37 Save your entries with **SAVE**.

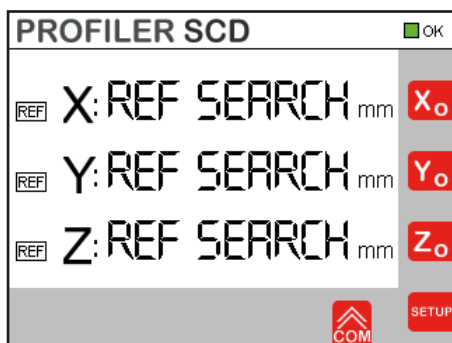38 Quit the menu *Encoder 1/2* with **ESC**.

39 Switch the PROFILER SCD off and on again.

Configuring the position correction is now completed.

In order to carry out corrected measurements, you only need to enable the position correction.

» *Chapter 6.5.6 Enabling the Position Correction, page 150*

## 6.5.4 Modifying Position Correction Values

### ► Modifying position correction values

- 1 Disable the *Corr* check box in the menu *Encoder 1/2* for the axis whose position correction values you want to modify.
- 2 Open the menu *Position Correction (PosCorr)* and switch to the menu *Position Correction 2/2* with **↓**.
- 3 Navigate through the saved segments with **▲** and **▼**.
- 4 To modify a target value, tap on the *Target Pos.* input field.

## NOTICE

### Incorrect position correction from excessive deviation between target value and actual value.

The *Target Pos.* and *Act. Pos.* values may deviate from each other by a maximum of 127  $\mu\text{m}$ . Check each value pair before saving it.

### Incorrect position correction from entering negative values.

Only positive values may be entered for *Target Pos.* and *Act. Pos.* Do not enter values with a negative sign.

- 
- 5 Enter the value of the measuring point approached on the measurement scale via the entry bar.
  - 6 Confirm your entry with .
  - 7 To modify an actual value, adopt the currently measured position value in the *Act. Pos.* input field with .
 

**Note:** If you would like to enter or modify the actual value manually, tap on the *Act. Pos.* input field. Enter or modify the actual value via the entry bar. Confirm your entry with .
  - 8 Save the entered target and actual value for the segment with .

The new position correction values are now saved.

- 9 Enable the *Corr* check box in the menu *Encoder 1/2* for the axis whose position correction values you have modified.

Modifying the position correction values is now completed. Before you can carry out corrected measurements, you only need to enable the position correction.

» *Chapter 6.5.6 Enabling the Position Correction, page 150*

## 6.5.5 Deleting Position Correction Values

### ► Deleting position correction values

- 1 To delete all segments which are saved for an axis along with the associated target and actual values, navigate with  to segment 0.
- 2 Delete all segments which were saved for the respective axis with **Erase**.
- 3 Confirm the query *Delete all position correction data of axis?* with **Yes**.

All segments which were saved for the respective axis, along with their associated target and actual values, are now deleted.

## 6.5.6 Enabling the Position Correction

**To enable the position correction, you must set/move to the machine zero** of your measuring stage/setup every time the PROFILER SCD is switched on.

When the machine zero is approached, the position correction status display of the X-, Y- and/or Z-axis changes to Green. The position correction is active. You can now carry out corrected measurements.

**Note:** If you do not set/move to the machine zero of an axis, the position display of the respective axis shows *REF SEARCH*; the *position correction status display* flashes red. In this case, the position correction is not enabled for the respective axis.

## 6.6 Enabling/Disabling the Standby Mode

In standby mode, the touchscreen of the PROFILER SCD is dimmed with the exception of the *general status display*. Power consumption is reduced. The measurement function is still active. The set machine zero and the connection to the PC are kept.

### ► Enable standby mode

- 1 Tap anywhere on the title bar of the position display.

The touchscreen turns black. Only the *general status display* changes to yellow.

### ► Disable standby mode

- 1 Tap on any part of the touchscreen.

The position display is shown again.

## 6.7 Configuring the PROFILER SCD

In the next chapters, you will learn how to configure the PROFILER SCD using different menus.

### 6.7.1 Setup

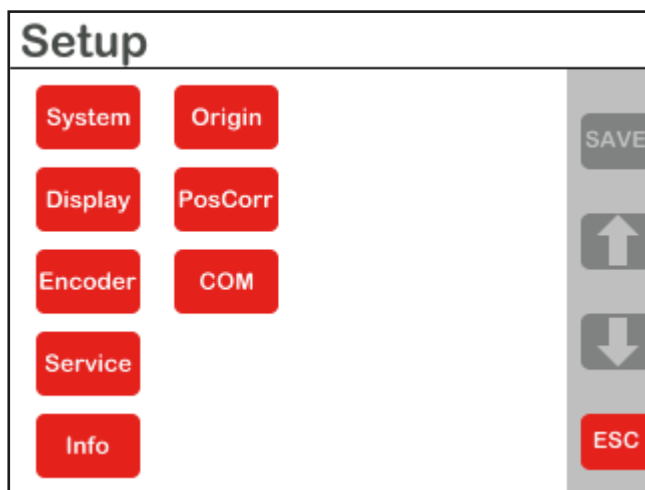
The menu *Setup* gives you access to all menus, in which you can define settings.

When password query is enabled, the menu *Setup* is protected with the password 1286. To make settings, enter the password and confirm your entry with .

You can disable the password query in the menu *System*.

» Chapter 6.7.2 *System*, page 153

#### **System**

configure operating properties

#### **Origin**

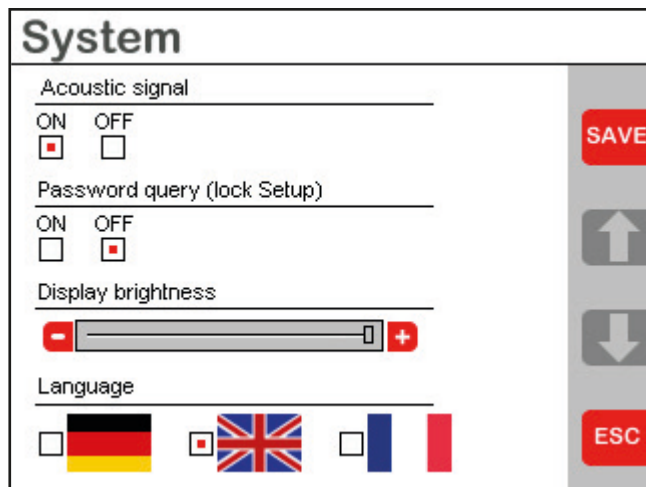
configure settings for zeroing the position value and for setting the machine zero

<b>Display</b>	configure display settings
<b>PosCorr</b>	configure position correction
<b>Encoder</b>	configure measuring system evaluation
<b>COM</b>	configure data transfer to a PC
<b>Service</b>	monitor signal quality of the measuring system, restore factory setting, perform reset
<b>Info</b>	view system information
<b>ESC</b>	return to position display

## 6.7.2 System

You can configure the basic operating properties of the PROFILER SCD in the menu *System*.

**Acoustic signal**      enable/disable acoustic signal

<b>Password query</b>	enable/disable password query
<b>Display brightness</b>	increase/decrease display brightness
<b>Language</b>	select user interface language
<b>SAVE</b>	save settings
<b>ESC</b>	return to menu <i>Setup</i>

When password query is enabled, the menu *Setup* is protected with the password 1286.

### 6.7.3 Origin

In the menus *Origin 1-3*, you can, separately for each axis (X, Y, Z),

- enable and disable the manual zeroing of the position values in the position display with **X<sub>o</sub>**, **Y<sub>o</sub>** and **Z<sub>o</sub>**,
- select the function for setting the machine zero, and
- enter the zero offset of the machine zero.

Distinguish between the following terms:

- **Position value:** Current position value, which is displayed for each axis in the position display, and can be zeroed with **X<sub>o</sub>**, **Y<sub>o</sub>** and **Z<sub>o</sub>**.
- **Machine zero:** Coordinate zero point of the respective axis. For each axis, the machine zero can be set manually or by a reference switch/reference mark.  
If you have set a function for setting the machine zero, the position display shows *REF SEARCH* for the respective axis as long as the machine zero is not set or approached.  
If you change the function for setting the machine zero, the current machine zero becomes lost and must be reset using the selected function.

Also distinguish between:

- **Reference switch:** In your measuring stage from Märzhäuser Wetzlar, integrated reference switches generate a control signal when they are switched, and thus the accurate reproducible machine zero for the position correction.
- **Reference mark:** Measuring systems from other manufacturers can be equipped with one or several reference marks, which deliver a separate control signal when they are approached. This control signal can also be used for setting the machine zero for the position correction.

## **NOTICE**

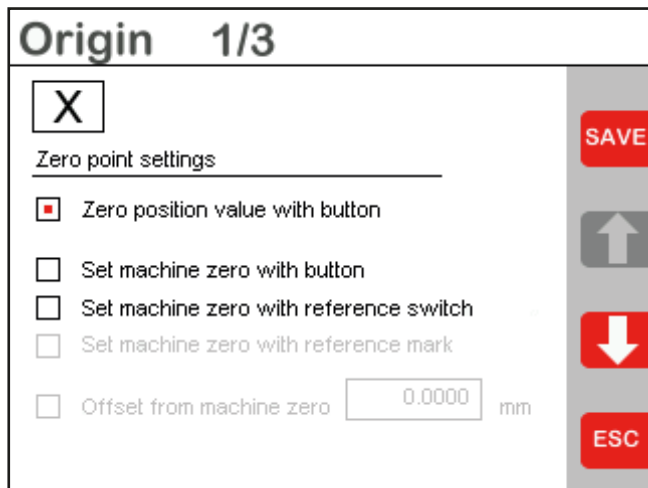
### **Incorrect position values from relocating the machine zero or changing the encoder type.**

Position correction values are valid as long as the location of the machine zero and the chosen encoder type are unchanged. If you, after entering position correction values,

- relocate the machine zero,
- approach another reference mark to set the machine zero,
- change the function for determining the machine zero or
- change the encoder type,

you must delete all previously stored position correction values and enter new position correction values.

---

Which functions are inactive in the menu *Origin* depends on the selected encoder type (for each axis) in the menu *Encoder*.

**Zero position value with button** enable/disable  $X_o$ ,  $Y_o$  and  $Z_o$  for the manual zeroing of position values

**Note:** The function *Zero position value with button* is only used for carrying out measurements independently of the machine zero.

» *Chapter 6.3 Performing Measurements, page 137*

**Set machine zero with button** set the machine zero with  $X_o$ ,  $Y_o$  and  $Z_o$

**Function description:** To set the machine zero, after enabling the function or after each time the PROFILER SCD is switched on, move your measuring stage to the desired position and manually set the machine zero one time with  $X_o$ ,  $Y_o$  and  $Z_o$ .

### Set machine zero with reference switch

$X_0$ ,  $Y_0$  and  $Z_0$  are then used again to zero the position values.

Reference signals are ignored. The position correction does not refer to the reference switches/reference marks, but to the manually set machine zero.

set the machine zero by approaching a reference switch

**Function description:** To set the machine zero using a reference switch, you must enable the reference signal evaluation for the respective axis in the menu *Encoder*.

» Chapter 6.7.6 Encoder, page 163

If the *REF* check box in the menu *Encoder* is enabled for an axis, the *status display reference signal* **REF** appears to the left of the respective position display.

To set the machine zero, approach the reference switch of the respective axis after enabling the function or after each start of the PROFILER SCD.

If a reference switch is approached, the *status display reference signal* **REF** to the left of the respective position display changes from white to green.

### Set machine zero with reference mark

set the machine zero of the 3rd axis by approaching the nearest reference mark

**Function description:** This function is available only for the 3rd measuring system, which is connected to the port *Encoder 3* of the PROFILER SCD (Z-axis, menu *Origin 3/3*).

To set the machine zero using a reference mark, you must enable the reference signal evaluation for the 3rd axis in the menu *Encoder*.

» *Chapter 6.7.6 Encoder, page 163*

If the *REF* check box in the menu *Encoder* is enabled for the 3rd axis, the *status display reference signal* **REF** appears to the left of the Z position display.

Move your 3rd axis to the desired position before you enable the function or switch on the PROFILER SCD. The machine zero is set with the first reference mark, which you approach after enabling the function or after switching on the PROFILER SCD.

If the reference mark is approached, the *status display reference signal* **REF** to the left of the position display Z changes from white to green.

The following reference marks are ignored.

**Note:** If your 3rd measuring system has several reference marks, be sure first to approach the reference mark, which you used for configuration of the position correction. Otherwise there can be no proper position correction.

**Offset from machine zero**

indicate the deviation between the machine zero and the zero point of the sample/workpiece

**Function description:** To exactly approach a determined zero point of your sample/workpiece after setting/approaching the machine zero, indicate the *Offset from machine zero* (deviation between machine zero and zero point of your sample/workpiece).

**Note:** The value can only be entered if you have selected a function for setting the machine zero.

**SAVE**

save settings

**Up/down arrow**

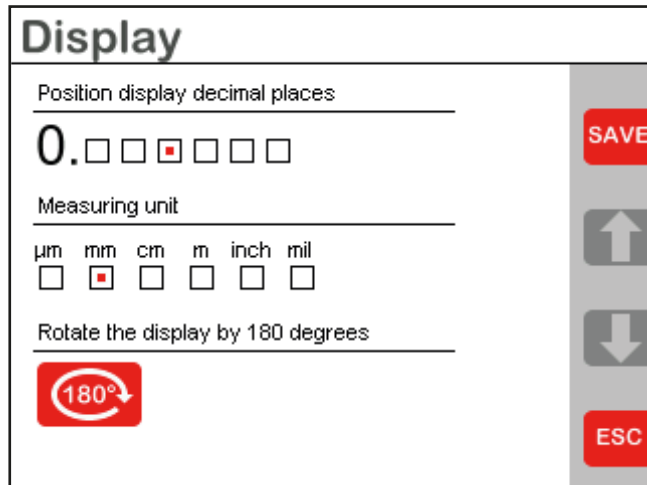
switch to the menu *Origin* of the other axes

**ESC**

return to the menu *Setup*

## 6.7.4 Display

You can configure the display settings of the PROFILER SCD in the menu *Display*.

**Position display decimal places**

define the number of displayed decimal places

**Measuring unit**

define measuring unit

**Rotate the display by 180 degrees**

rotate touchscreen display by 180° (e.g. for wall installation)

**SAVE**

save settings

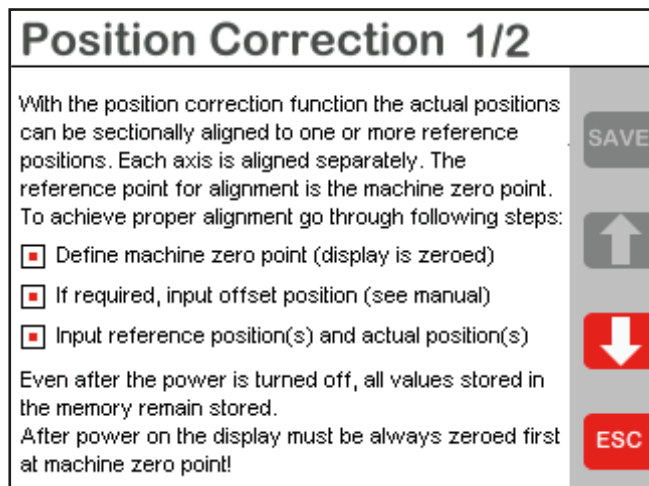
**ESC**

return to the menu *Setup*

## 6.7.5 PosCorr (Position Correction)

The menu *Position Correction 1/2* describes the procedure for configuring the position correction.

**Down arrow**            switch to the menu  
*Position Correction 2/2*

**ESC**                    return to the menu *Setup*

In the menu *Position Correction 2/2*, you can configure the position correction of the PROFILER SCD.

Position Correction 2/2		
<b>X</b>	<b>Y</b>	<b>Z</b>
Target Pos. 13.000	Target Pos. 0.000	Target Pos. 0.000
Act. Pos. <b>L</b> 12.475	Act. Pos. <b>L</b> 0.000	Act. Pos. <b>L</b> 0.000
Segment 1	Segment 0	Segment 0
<b>Enter</b>	<b>Erase</b>	<b>Erase</b>

- Target Pos.** enter the target value of the reference position
- Act. Pos.** enter the actual value of the reference position
- L** display live the current position value in the *Act. Pos.* field
- Segment** segment number, to which the displayed target and actual values are assigned
- Enter** save the displayed/entered target and actual values from the fields *Target Pos.* and *Act. Pos.* for the displayed segment
- Erase** delete target and actual values of all segments of an axis
- Note:** **Erase** is only visible when segment 0 is selected.

<b>Up arrow</b>	switch to the menu <i>Position Correction 1/2</i>
<b>ESC</b>	return to the menu <i>Setup</i>

Here you can find information on configuring and enabling the position correction:

» *Chapter 6.5 Using the Position Correction, page 140*

## 6.7.6 Encoder

In the menu *Encoder 1/2*, you can select different evaluation options for the connected measuring systems.

### **NOTICE**

#### **Incorrect measurement from wrong configuration.**

When configuring the PROFILER SCD, you must indicate the correct encoder type for the 3rd measuring system, if available.

#### **Incorrect position values from relocating the machine zero or changing the encoder type.**

Position correction values are valid as long as the location of the machine zero and the chosen encoder type are unchanged. If you, after entering position correction values,

- relocate the machine zero,
- approach another reference mark to set the machine zero,
- change the function for determining the machine zero or
- change the encoder type,

you must delete all previously stored position correction values and enter new position correction values.

Encoder 1/2		
<b>X</b>	<b>Y</b>	<b>Z</b>
<input checked="" type="checkbox"/> ON/OFF	<input checked="" type="checkbox"/> ON/OFF	<input checked="" type="checkbox"/> ON/OFF
<input checked="" type="checkbox"/> Count +/-	<input checked="" type="checkbox"/> Count +/-	<input checked="" type="checkbox"/> Count +/-
<u>Encoder type</u>	<u>Encoder type</u>	<u>Encoder type</u>
MR <input checked="" type="checkbox"/> TTL <input type="checkbox"/> 1Vss <input type="checkbox"/>	MR <input checked="" type="checkbox"/> TTL <input type="checkbox"/> 1Vss <input type="checkbox"/>	MR <input type="checkbox"/> TTL <input checked="" type="checkbox"/> 1Vss <input type="checkbox"/>
REF <input type="checkbox"/> Corr <input type="checkbox"/>	REF <input type="checkbox"/> Corr <input type="checkbox"/>	REF <input type="checkbox"/> Corr <input type="checkbox"/>
Period [mm] 0.500000	Period [mm] 0.500000	Counting inc. 0.00025
<input type="button" value="SAVE"/> <input type="button" value="↑"/> <input type="button" value="↓"/> <input type="button" value="ESC"/>		

**ON/OFF**

enable/disable measurement function

**Note:** When the measurement function is disabled, the respective axis is hidden in the position display.

**Count +/-**

select counting direction

**Encoder type**

select encoder type  
(MR/analog 5Vss, 1Vss, TTL (RS-422))

**Note:** The selection of the encoder types 1Vss and TTL (RS-422) is only possible for the 3rd measuring system.

**REF**

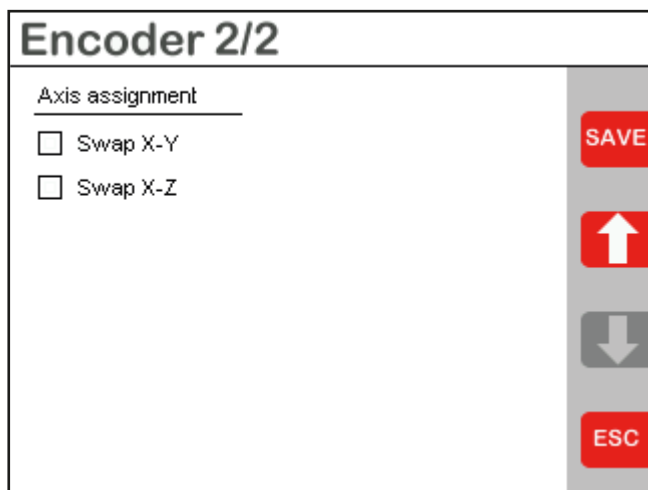
enable/disable reference signal evaluation

**Corr**

enable/disable position correction


<b>Period</b>	<i>MR/analog 5Vss</i> : 0.5 mm pre-set (not adjustable) <i>1Vss</i> : enter period length of the input signal in mm <i>TTL (RS-422)</i> : enter counting step according to quadruple evaluation in mm
<b>SAVE</b>	save settings
<b>Down arrow</b>	switch to the menu <i>Encoder 2/2</i>
<b>ESC</b>	return to the menu <i>Setup</i>

In the menu *Encoder 2/2*, you can swap the axis assignment of the connected measuring systems.

### Axis assignment

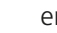enable/disable swap of the position displays for the X- and Y-axis or the X- and Z-axis

**Note:** When *Swap X-Y* or *Swap X-Z* is enabled, arrows  show which axes are switched in the menu *Encoder 1/2*. The setting *Swap X-Y* or *Swap X-Z* does not affect the assignment of reference switches/marks and position correction values to the related axes. Only the display of the position values on the PROFILER SCD is switched.

### SAVE

save settings

### Up arrow

switch to the menu *Encoder 1/2*

### ESC

return to the menu *Setup*

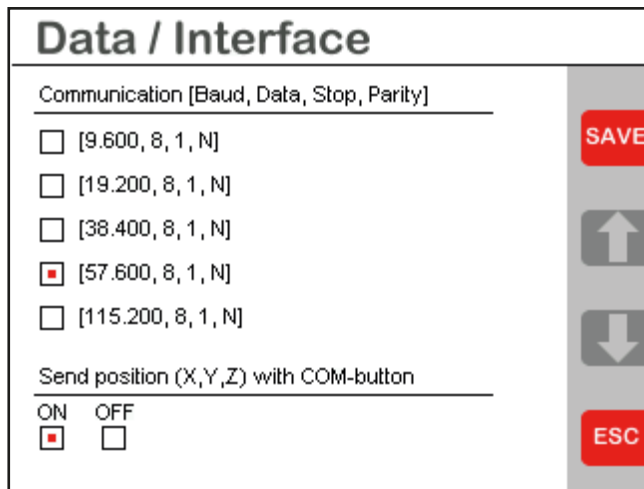
### 6.7.7 COM (Data / Interface)

In the menu *COM (Data / Interface)* you can configure the serial PC interface of the PROFILER SCD and enable/disable the button .

#### NOTICE

#### Faulty data transmission and measurement from incorrect configuration.

When the PROFILER SCD and PC are connected to each other over the RS-232 interface, you must enter the same baud rate selected on your PC when configuring the PROFILER SCD.

#### Communication

select baud rate

**Note:** The baud rate selection is only relevant for an RS-232 connection to your PC. With a USB connection, the baud rate selection has no effect.

<b>Send position (X, Y, Z) with COM-button</b>	enable/disable button 
<b>SAVE</b>	save settings
<b>ESC</b>	return to the menu <i>Setup</i>

## 6.7.8 Service

In the menu *Service 1/2*, you can see the current status of the connected measuring systems.

	Encoder X	Encoder Y	Encoder Z	
Position:	12.465	-5.148	14.740	SAVE
Hw. Count:	158	-64	184	↑
Amplitude:	84%	82%	78%	↓
Sin:	-111	144	75	ESC
Cos:	-404	-347	-402	
Offset Sin:	84	42	24	
Offset Cos:	-12	-3	-13	
Cos Amplify:	1.004	1.008	1.002	

**Down arrow** switch to the menu *Service 2/2*

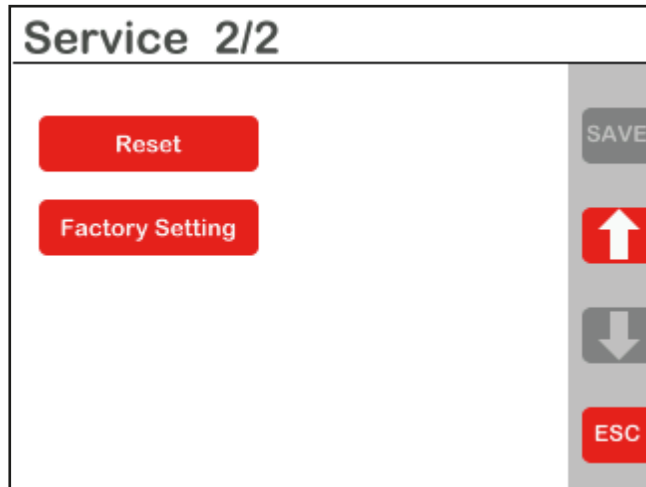
**ESC** return to the menu *Setup*

The values displayed in the menu *Service 1/2* are mainly for service purposes. You could be asked for the values which are displayed here when you have contact with our Customer Service Center in case of any problems.

The menu *Service 2/2* gives you the opportunity to perform a reset and to restore the factory setting of the PROFILER SCD.

### Reset

restart PROFILER SCD

**Note:** The saved settings are kept. After resetting, you only need to set/ approach the machine zero of your measuring stage, if the position correction is enabled.

### Factory Setting

reset PROFILER SCD to factory setting

**Note:** Saved position correction values remain kept when restoring the factory setting.

### Up arrow

switch to the menu *Service 1/2*

### ESC

return to the menu *Setup*

## 6.7.9 Information

In the menu *Information*, you can find information on the firm-ware version and type of the PROFILER as well as our contact data.

**ESC**

return to the menu *Setup*

## 7 Operation on a PC

### 7.1 GetPos Software

Using the *GetPos* software, you can send position values to your PC, and have the option of saving them in an ASCII text file (.txt) or an Excel-compatible file (.csv) for further processing.

To send position values from the PROFILER SCD to your PC, and to process them with *GetPos*, the PROFILER SCD and your PC must be correctly connected and configured.

» *Chapter 5.4 Establishing the Connection to a PC, page 126*

#### 7.1.1 Installation and Configuration

##### ► Installing *GetPos*

- 1 Create a new folder on your hard drive, to which you would like to install *GetPos*. Name the folder, e.g. *GetPos*.
- 2 Insert the installation CD in the CD drive.
- 3 Open the `..\Software\GetPos` folder.
- 4 Copy the file *GetPos.exe* to the folder on your hard drive which you created in step 1.

*GetPos* is now installed on your PC.

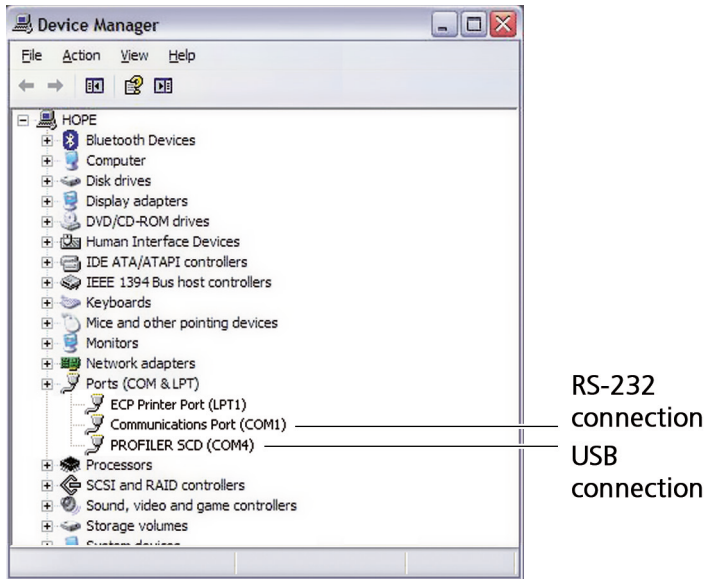
When using *GetPos* for the first time, it generates the configuration file *GetPos.cfg*, in which the settings adopted by you are saved. The configuration file is saved in the folder, in which the file *GetPos.exe* is stored.

##### ► Identifying the COM port

To configure *GetPos*, you need to identify the COM port, which is used by the PROFILER SCD.

- 1 Open the Device Manager:
  - **Windows 7:** *Start* → *Control Panel* → *System* → *Device Manager*
  - **Windows 8:** right-click in the left, lower corner of the desktop/start screen → *Device Manager*
  - **Windows 10/11:** right-click on the Windows symbol in the left, lower corner of the desktop/start screen → *Device Manager*
- 2 Expand the column *Ports (COM and LPT)* by double-clicking.

The COM port used is listed in brackets behind the entry *PROFILER SCD* (with a USB connection) or behind the entry *Communications Port* (with an RS-232 connection).


**RS-232 connection:** If your PC has more than one serial port, refer to the operating manual of your PC for information about the assignment of COM ports to serial ports.

**USB connection:** The number of the virtual COM port is kept regardless of which USB port of your PC you are using.

## ► Configuring GetPos

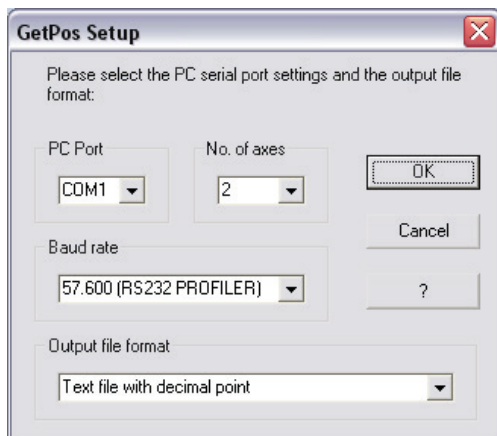- 1 Start GetPos by double-clicking on the file *GetPos.exe*.

When starting GetPos for the first time, the dialog window *GetPos Setup* appears. Here, you can make all required settings, in order to establish the connection between PROFILER SCD and GetPos.

If the settings were already made, the dialog window *GetPos Setup* no longer appears when starting GetPos. You can, however, adjust all settings any time over the menu *Setup* in the menu bar.

» *Chapter 7.1.3 Menu Functions, page 176*

## NOTICE

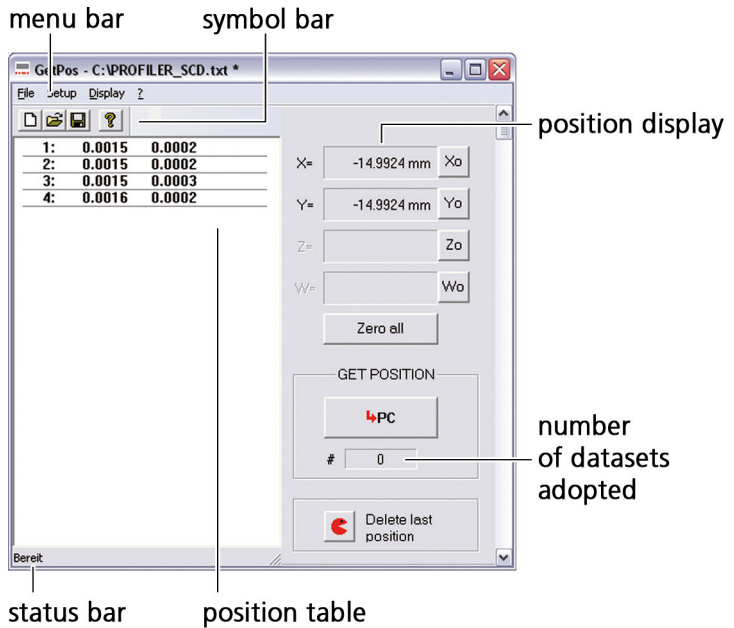
### Faulty data transmission and measurement from incorrect configuration.

When configuring GetPos, you must enter the correct port, the correct number of measuring systems (axes) and, when connected over the RS-232 interface, the baud rate selected on the PROFILER SCD.

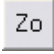
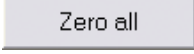
- 2 Select the *PC Port* over which the PROFILER SCD is connected to your PC.
  - 3 Select the number of measuring systems (*No. of axes*), which are connected to the PROFILER SCD and are enabled in the menu *Encoder*.
    - » *Chapter 6.7.6 Encoder, page 163*
  - 4 Select the same *Baud rate*, which is selected in the PROFILER SCD (factory setting: 57,600 baud).
    - » *Chapter 6.7.7 COM (Data / Interface), page 167*
- Note:** The baud rate selection is only relevant for an RS-232 connection to your PC. With a USB connection, the baud rate selection has no effect.
- 5 Select the *Output file format* which should be used when saving your data.
  - 6 Confirm your entries by clicking on the button *OK*.

GetPos is now configured for communication with the PROFILER SCD.

## 7.1.2 User Interface

Button	Function
	create new position table
	open saved position table
	save current position table
	display version number of GetPos
	zero position value of X

Button	Function
	zero position value of Y
	zero position value of Z
	no function <sup>1</sup>
	zero position values of all axes
	adopt current position values in position table
	delete last adopted position value

<sup>1</sup> function only available for predecessor devices with 4 measuring channels

### 7.1.3 Menu Functions

Menu	Entry	Function
File	New	create new position table
	Open...	open saved position table
	Save	save current position table
	Save as...	save current position table with a new filename

Menu	Entry	Function
	Last file	open last saved position table
	Quit	end GetPos
Setup	COM port	select port over which the PROFILER SCD is connected to your PC
	Baud rate	select baud rate <b>Note:</b> The baud rate selection is only relevant for an RS-232 connection to your PC. With a USB connection, the baud rate selection has no effect.
	File format	select file format in which the position table is saved
	No. of axes	select number of measuring systems, which are connected to the PROFILER SCD
	Language	select language (German/English)
Display	Tool bar	show/hide symbol list
	State bar	show/hide status bar
?	About Get-Pos...	display version number of GetPos

## 7.1.4 Key Layout on a PC

Keys/ key combinations	Function
X Y Z	zero position value of X, Y, Z
0	zero position values of all axes
ENTER	send current position values to PC
CTRL + Backspace	delete last adopted position values

## 7.1.5 Sending Position Values to the PC

You can send position values to GetPos in different ways:

- press button  on PROFILER SCD
- press button  in GetPos
- press button *Enter* on your PC keyboard
- press right pedal of the foot switch (optional accessory)

When you have sent the position values in one of the ways described, the position values appear in the GetPos position table as a new data set.

## 7.1.6 Saving Position Tables

You can save position tables in the following file formats:

- text file with decimal separators (ASCII)
- text file with comma separators (ASCII)
- Excel-compatible file (.csv)

### ► Changing the file format

When starting GetPos for the first time, you determine different settings, including the file format, in which you will save your position tables. These settings can, however, be adjusted any time over the menu *Setup* in the menu bar.

- 1 In the menu *Setup* in the menu bar, select the preferred file format in the sub menu *File format*.

The position table is saved in the selected file format the next time it is saved.

### ► Saving a position table

- 1 In the menu *File* in the menu bar, select the entry *Save* or click on  in the tool bar.
- 2 Follow the on-screen instructions.

### ► Saving a position table with a new filename

- 1 In the menu *File* in the menu bar, select the entry *Save as...*
- 2 Follow the on-screen instructions.

### ► Open a position table

GetPos automatically creates a new position table with each program start.

- 1 To open a saved position table, click on the button  in the tool bar or select the entry *Open...* in the menu *File* in the menu bar.
- 2 Follow the on-screen instructions.

As soon as the position table is visible, you can add additional position values and save the file under a new name or in a different file format.

## 7.2 Instruction Set

Using instruction sets, you can read position values and configure all settings of the PRFOILER SCD over the RS-232 or USB interface.

The following instruction sets are available:

- **Instruction Set PROFILER**  
The original instruction set of the PROFILER product family provides comprehensive options for the programming of the PROFILER SCD. The use is recommended.
- **Instruction Set SCD<sup>plus</sup>** (applied for PROFILER SCD)  
The instruction set is compatible with the instruction set of the predecessor devices *SCD5*, *SCD6* and *SCD<sup>plus</sup>* (with few exceptions) and is suitable for programming the PROFILER SCD, if you want to use or develop further an already existing software.

You can find the instruction sets available in PDF format on the installation CD:

- » *Instruction Set PROFILER*
- » *Instruction Set SCD<sup>plus</sup>*

## 8 Cleaning

### WARNING

#### **Danger to life and risk of equipment damage from improper cleaning.**

Disconnect the power supply from the mains before cleaning the PROFILER SCD or the power supply. Ensure that no liquids enter the interior of the PROFILER SCD or the power supply. Protect ports and plugs from liquids. Never touch the power supply or the socket with wet hands.

#### **Danger to life and risk of equipment damage from spreading hazardous substances.**

In case of possible contamination with corrosive, infectious, toxic, radioactive or other substances which are harmful to health, have the PROFILER SCD professionally disinfected/decontaminated.

---

#### ► **Cleaning the PROFILER SCD**

- 1 Disconnect the PROFILER SCD from the mains.
- 2 Clean the PROFILER SCD with a lint-free cleaning cloth. With greater contamination, slightly dampen the cleaning cloth with water or a mild solvent-free disinfectant/cleaning agent.
- 3 Allow the PROFILER SCD to dry.
- 4 Connect the PROFILER SCD to the mains.

The PROFILER SCD can now be used again.

## **9 Maintenance**

The PROFILER SCD is maintenance-free.

## 10 Troubleshooting

### WARNING

#### **Possible electric shock and equipment damage from connecting foreign or damaged power supplies.**

Only use the delivered 24 V/30 W power supply. Immediately replace the power supply if there is damage on the housing or cable of the power supply. Only use replacement components from Märzhäuser Sensotech.

#### **Possible electric shock and equipment damage from not permitted repair attempts.**

Never open the housing of the PROFILER SCD or the accompanying power supply. Only have repair work carried out by our Customer Service Center. In case of violation, all warranty and liability claims will become void.

---

The following overview can help you in removing faults and malfunctions of the PROFILER SCD.

If a fault or malfunction cannot be remedied using the overview, contact our Customer Service Center.

» *Chapter 11 Service, page 188*

<b>Fault</b>	<b>Possible causes</b>
PROFILER SCD doesn't turn on	no power supply
Position display shows <i>NO SENSOR</i> for an axis	incorrect cabling
	measuring channel is not needed
Position display shows <i>REF SEARCH</i> for an axis	machine zero not set/approached
Position values unusual or demonstrably incorrect	incorrect information about the encoder type or signal period/ counting step
	faulty position correction values (possible when you have entered or modified position correction values yourself)
	connection of an incompatible measuring system

## Remedy

Check whether the external power supply is properly connected to the port 24V/DC on the PROFILER SCD and the socket.

- 1 Disconnect the power supply of the PROFILER SCD from the mains.
- 2 Check whether the affected measuring system is properly connected to the PROFILER SCD over the correct port.
- 3 Connect the power supply of the PROFILER SCD to the mains.
- 4 Check the position display of the affected axis.

Disable the position displays of unused measuring channels.

» *Chapter 6.7.6 Encoder, page 163*

- If function *Set machine zero with button* is selected: After switching on the PROFILER SCD, manually set the machine zero with **X<sub>0</sub>**, **Y<sub>0</sub>** or **Z<sub>0</sub>**.
- If function *Set machine zero with reference switch* or function *Set machine zero with reference mark* is selected: Approach the reference switches/reference marks of all axes after switching on the PROFILER SCD.

» *Chapter 6.7.3 Origin, page 154*

Check the settings for the encoder type and the signal period/counting step of the affected measuring system.

» *Chapter 6.7.6 Encoder, page 163*

- 1 Check the position correction values of the affected measuring system.
- 2 Correct faulty position correction values or, if necessary, delete all position correction values and configure the position correction once more.

» *Chapter 6.5.3 Configuring the Position Correction Yourself, page 141*

Check the specifications of the affected measuring system.

» *Chapter 4.1 Intended Use, page 116*

<b>Fault</b>	<b>Possible causes</b>
Position correction does not work	machine zero not set/ approached (position display shows <i>REF SEARCH</i> , <i>status display position correction</i> of the respective axis flashes red)
	position correction for the affected axis is disabled

## Remedy

---

- If function *Set machine zero with button* is selected: After switching on the PROFILER SCD, manually set the machine zero with **X<sub>0</sub>**, **Y<sub>0</sub>** or **Z<sub>0</sub>**.
  - If function *Set machine zero with reference switch* or function *Set machine zero with reference mark* is selected: Approach the reference switches/reference marks of all axes after switching on the PROFILER SCD.
    - » *Chapter 6.7.3 Origin, page 154*
- 

Enable the position correction for the affected axis.

- » *Chapter 6.7.6 Encoder, page 163*
-

## 11 Service

The staff at our Customer Service Center are happily at your disposal for support queries, complaints and suggestions about our products and services.

Overview of the services from our Customer Service Center:

- advice and support
- support for customer-specific application questions
- support for putting equipment into service
- inspection and repair of equipment
- error analysis (also online via web conference)
- delivery of replacements and replacement components
- user training (also online via web conference)

You can reach our Customer Service Center Mondays to Fridays from 08:00 am to 04:00 pm CET.

Tel.: **+49 6441 9116-36**  
 Fax: **+49 6441 9116-40**  
 E-mail: **[service@marzhauser.com](mailto:service@marzhauser.com)**

So that we can help you quickly when you need service, please keep handy the following information about your PROFILER SCD:

Description	Information on the identification plate
Type	
Article-No.	
Serial-No.	

## 11.1 Replacement Components

Should you need to replace a component of the PROFILER SCD, please contact our Customer Service Center.

» *Chapter 11 Service, page 188*

<b>Replacement components</b>	<b>Order no.</b>
MR encoder cable (2 channels, 2 m)	00-99-900-0859
RS-232 interface cable	00-76-232-0801
USB interface cable	00-76-100-2802
External 24 V/30 W power supply incl. 4 adapters (EU/GB/USA/AUS)	00-76-203-4801
Operating manual	00-76-900-4830-TD
Installation CD incl. driver, oper- ating manual, instruction set, software <i>GetPos</i>	00-76-900-4830-CD

## 12 Taking out of Service

### ► Taking the PROFILER SCD out of service

- 1 Switch off the PROFILER SCD (switch *ON-OFF* to position *0*).

### **NOTICE**

#### **Possible equipment damage from putting into service / taking out of service with existing power connection.**

Disconnect the PROFILER SCD from the mains before connecting/disconnecting measuring systems, a PC or other peripherals to/from the PROFILER SCD.

- 2 Disconnect the power supply of the PROFILER SCD from the mains.
- 3 Disconnect all connected cables from the PROFILER SCD.
- 4 Clean the PROFILER SCD.  
» *Chapter 8 Cleaning, page 181*

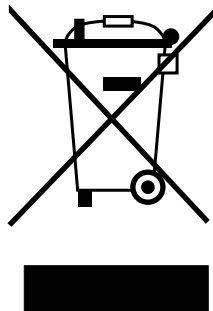
The PROFILER SCD can now be disposed of.

## 13 Disposal

### NOTICE

#### Possible punishable act through improper disposal.

In the EU, the disposal of electrical and electronic equipment with your household waste is illegal. Do not dispose of the PROFILER SCD and accessories with the household waste. If you dispose of the PROFILER SCD outside the EU, note the legal regulations.

The adjacent symbol is attached to the PROFILER SCD. This symbol means that the PROFILER SCD is classified based on the existing regulations as electrical and electronic equipment, and falls under EU directive 2012/19/EU (WEEE). The PROFILER SCD may not, therefore, be disposed of with household waste.

Märzhäuser Sensotech will properly dispose of the PROFILER SCD.

#### ► Disposing of the PROFILER SCD

- 1 Take the PROFILER SCD out of service.
  - » *Chapter 12 Taking out of Service, page 190*
- 2 Properly pack the PROFILER SCD and accessories.
- 3 Send the PROFILER SCD for disposal to:

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### 14.1 General Data

#### Product features

Measuring channels	3
Position resolution	up to 1 nm (depends on measuring system used)
Interpolation accuracy	0.1 % of the signal period (MR/analog 5Vss, 1Vss)
Signal inputs	2 × MR/analog 5Vss 1 × configurable: MR/analog 5Vss, 1Vss, TTL (RS-422)
Specification of the signal inputs	MR/analog 5Vss, 1Vss: interpolation up to factor 12,868 TTL (RS-422): up to 25 MHz
Max. travel speed (X, Y)	2.5 m/s
Max. input frequency (Z)	25 MHz
USB interface	USB 2.0 Full Speed (12 Mbit/s)
RS-232 interface	signals RXD/TXD, up to 115 kbaud
Feature interface	optional functions (port for foot switch, reference switch input for 3rd measuring system)
Programming	ASCII

### Product features

Touchscreen	320 × 240 pixels, resistive
Housing	aluminium, plastics
Protection class	II
Protection type	IP20
Supply voltage	8...28 V DC, typ. 110 mA at 24 V (without connected peripherals)
Power consumption	typ. 4 W (operation) and 2 W (standby mode) with 2 connected measuring systems
Dimensions	210 × 190 × 65 mm (L × W × H)
Weight	approx. 0.9 kg (excluding cables)

On request, your PROFILER SCD can also be upgraded for the connection of an incremental measuring system of type *TTL (non-differential)* on the port *Encoder 3*. When needed, please contact our Customer Service Center:

» *Chapter 11 Service, page 188*

### System requirements (when connected to a PC)

Interface	USB or RS-232 interface
Operating system	Windows 7 (32/64 bit), Windows 8 (32/64 bit), Windows 10 (32/64 bit) or Windows 11 (64 bit)

### Security functions

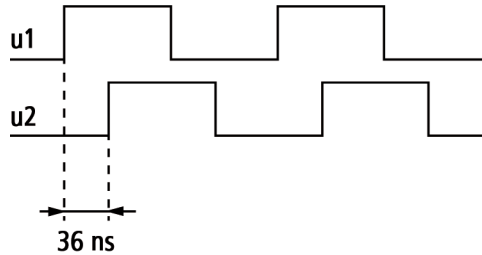
Processor monitoring	reset in case of undervoltage, logic voltage (3.3 V)
Voltage monitoring	measurement of internal voltages (3.3 V, 5 V, 24 V/DC)
Surge/short-circuit protection (external voltages)	VCC5 + VCCEXT, shutdown is automatically reset after the end of the overload condition

## 14.2 Encoder Interface

Parameter	MR/ana- log 5Vpp	1Vss	TTL (RS-422)
Position signal	1.5...5 Vss <sup>1</sup>	0.6...1.2 Vss <sup>1</sup>	standard <sup>1</sup>
Reference signal evaluation	–	yes	yes
Error signal evaluation	on request	on request	on request
Max. input frequency	25 MHz	25 MHz	25 MHz <sup>2</sup>
Max. interpolation factor	12,868 <sup>3</sup>	12,868 <sup>3</sup>	4 <sup>4</sup>
Input resistor	120 Ω	120 Ω	120 Ω
Supply voltage measuring system	5 V <sup>5</sup>	5 V <sup>5</sup>	5 V <sup>5</sup>

<sup>1</sup> differential transmission

<sup>2</sup> pulse interval min. 36 ns

<sup>3</sup> arcus tangent interpolation

<sup>4</sup> quadrature counter

<sup>5</sup> max. 500 mA per measuring system, max. 800 mA total

## 14.3 Ports

On all ports, the cable shield is connected to the housing of the PROFILER SCD.

### 14.3.1 Encoder 1: HD D-Sub 15-Pin Female Connector

Pin	Signal	Function
1, 2	VCC5	power supply, 5 V <sup>1</sup>
3	+U1X1	+sine, X-axis <sup>2</sup>
4	-U1X1	-sine, X-axis <sup>2</sup>
5	+U2X1	+cosine, X-axis <sup>2</sup>
6	-U2X1	-cosine, X-axis <sup>2</sup>
7	REF2	reference switch input Y-axis <sup>3</sup>

Pin	Signal	Function
8	REF1	reference switch input X-axis <sup>3</sup>
11	+U1X2	+sine, Y-axis <sup>2</sup>
12	-U1X2	-sine, Y-axis <sup>2</sup>
13	+U2X2	+cosine, Y-axis <sup>2</sup>
14	-U2X2	-cosine, Y-axis <sup>2</sup>
9, 10, 15	GND	reference potential

<sup>1</sup> max. 500 mA, total load of all VCC5 ports max. 800 mA

<sup>2</sup> output signal: 5V<sub>ss</sub>, differential transmission, R<sub>L</sub> = 120 Ω

<sup>3</sup> TTL input, pull-up resistor 1 kΩ → VCC5, low pass 1 kΩ/10 nF

### 14.3.2 Encoder 3: D-Sub 9-Pin Female Connector

#### Pin assignment for MR/analog 5Vpp measuring systems

Pin	Signal	Function
1	-U1	-sine <sup>1</sup>
2	GND	reference potential
3	-U2	-cosine <sup>1</sup>
4	ERR	error <sup>2</sup>
5	-	-
6	+U1	+sine <sup>1</sup>
7	VCC5	power supply, 5 V <sup>3</sup>

Pin	Signal	Function
8	+U2	+cosine <sup>1</sup>
9	–	–

<sup>1</sup> output signal: 5V<sub>ss</sub>, differential transmission, R<sub>L</sub> = 120 Ω

<sup>2</sup> TTL input, pull-down resistor 1 kΩ → GND,  
low pass 1 kΩ/10 nF

<sup>3</sup> max. 500 mA, total load of all VCC5 ports max. 800 mA

### Pin assignment for 1V<sub>ss</sub> measuring systems

Pin	Signal	Function
1	–U1	–sine <sup>1</sup>
2	GND	reference potential
3	–U2	–cosine <sup>1</sup>
4	ERR	error <sup>2</sup>
5	–U0	–reference mark measuring system <sup>4</sup>
6	+U1	+sine <sup>1</sup>
7	VCC5	power supply, 5 V <sup>3</sup>
8	+U2	+cosine <sup>1</sup>
9	+U0	+reference mark measuring system <sup>4</sup>

<sup>1</sup> output signal: 1V<sub>ss</sub>, differential transmission, R<sub>L</sub> = 120 Ω

<sup>2</sup> TTL input, pull-down resistor 1 kΩ → GND,  
low pass 1 kΩ/10 nF

<sup>3</sup> max. 500 mA, total load of all VCC5 ports max. 800 mA

<sup>4</sup> differential transmission, R<sub>L</sub> = 120 Ω

**Pin assignment for TTL (RS-422) measuring systems**

Pin	Signal	Function
1	-U1	negative U1 <sup>1</sup>
2	GND	reference potential
3	-U2	negative U2 <sup>1</sup>
4	ERR	error <sup>2</sup>
5	-U0	-reference mark measuring system <sup>1</sup>
6	+U1	positive U1 <sup>1</sup>
7	VCC5	power supply, 5 V <sup>3</sup>
8	+U2	positive U2 <sup>1</sup>
9	+U0	+reference mark measuring system <sup>1</sup>

<sup>1</sup> output signal: TTL, differential transmission,  $R_L = 120 \Omega$

<sup>2</sup> TTL input, pull-down resistor  $1 \text{ k}\Omega \rightarrow \text{GND}$ ,  
low pass  $1 \text{ k}\Omega/10 \text{ nF}$

<sup>3</sup> max. 500 mA, total load of all VCC5 ports max. 800 mA

### 14.3.3 RS-232: D-Sub 9-Pin Male Connector

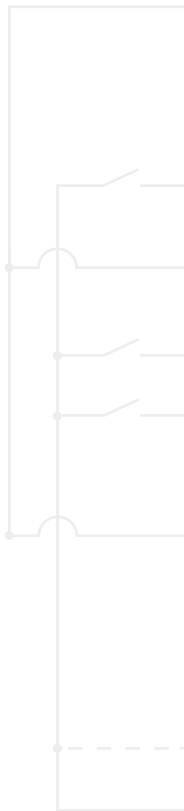
Pin	Signal	Function
1	–	–
2	RXD	reception line
3	TXD	transmission line
4	–	–
5	GND	reference potential
6	–	–
7	–	–
8	–	–
9	–	reserved for system-internal functions

### 14.3.4 USB Port: 4-Pin Female Connector, Type B

Pin	Signal	Function
1	VCC <sup>1</sup>	power supply (+5 V, max. 500 mA)
2	USBDM	–Data
3	USBDP	+Data
4	GND	reference potential

<sup>1</sup> internal 100 nF → GND

### 14.3.5 Feature: HD D-Sub 15-Pin Male Connector

Pin	Signal	Function
1	–	–
2	–	–
3	GND	reference potential
4	–	–
5	–	–
6	+IN1	+reference switch, 0-24 V <sup>1</sup>
7	-IN1	-reference switch, 0-24 V <sup>1</sup>
8	IN2	foot switch 1, Clear, 0-24 V <sup>1</sup>
9	IN3	foot switch 2, Send, 0-24 V <sup>1</sup>
10	–	–
11	–	reference potential IN2, IN3
12	NC	not connected
13	–	reserved for system-internal functions
14	VCC5	output voltage 5 V <sup>2</sup>
15	VCCEXT	output voltage 24 V ± 10% <sup>3</sup>

<sup>1</sup> logic level: 0–0.8 V = low; 3.95–28 V = high; secured against reverse polarity, optically isolated

<sup>2</sup> max. 500 mA, total load of all VCC5 ports max. 800 mA

<sup>3</sup> max. 0.4 A

### 14.3.6 24V/DC: 2.1 mm DC Male Connector

Pin	Signal	Function
Pin, inside	+24 V	voltage range +8... +28 V, power consumption at +24 V = typ. 110 mA, no load, internal fuse (not replaceable) = 2 A <b>Note:</b> Only use the delivered 24 V/30 W power supply.
Housing, external	GND	shield connection (connected to the housing)

### 14.3.7 Shield: Insert Nut M3

On the insert nut M3 (*Shield*), you can set the PROFILER SCD to a specific potential (e.g. protective earth).

The port is not suitable for implementing a standards-compliant PE connection.

# 14.4 External 24 V/30 W Power Supply (Phihong PSAA30R-240)

## 30W Interchangeable AC Clip Adapter

PSAA30R-XXX-R Series

● 12V/15V/24V/56V ● DoE VI ● Interchangeable AC Clip ● Cable ● Black

The 30W interchangeable AC clip adaptor series is the latest offering from a world leader in charging adapter technology and manufacturing - Phihong Technology. Phihong 30W interchangeable AC clip adaptor series are designed to work with most 12V/15V/24V/56V powered devices and conform to most of the industry standards and meet DoE level VI energy efficiency requirements. We ensure our products meet or exceed all the industry performance and safety requirements.

### Features :

<b>Efficiency</b>	DoE Level VI (86.95% avg)
<b>Power Saving</b>	<0.1W @230Vac/50Hz
<b>Protection</b>	SCP, OVP, OCP
<b>Turn On Delay Time</b>	<3s @ 115Vac, max load
<b>Hold-up Time</b>	>10ms @ 115Vac, max load
<b>Rise Time</b>	<100ms @ 115Vac, max load
<b>Isolation</b>	Primary to Secondary 3000Vac, 10mA for 1min.
<b>Insulation</b>	Primary to Secondary >7M ohm 500Vdc
<b>DC Output Cable (12V)</b>	1500±50mm, UL2468, AWG#18
<b>(15V, 24V, 56V)</b>	1500±50mm, UL2468, AWG#24
<b>DC Output Connector</b>	5.5x2.1x10mm, center positive, straight barrel
<b>Dimension</b>	82.7*55*39.1 (mm)
<b>Weight</b>	180 (g)

### Electrical Specifications :

<b>Input Voltage</b>	100~240Vac (90~264Vac min/max)
<b>Frequency</b>	50/60Hz Nom (47~63Hz min/max)
<b>Input Current</b>	0.8A (RMS) @ 120Vac max load 0.5A (RMS) @ 240Vac max load
<b>Output Voltage</b>	12V / 15V / 24V / 56V
<b>Output Current</b>	2.5A / 2A / 1.25A / 0.54A
<b>Maximum Output Power</b>	30W continuous
<b>Output Ripple &amp; Noise</b>	≤ 120mV (pk-pk) for 12V output
<b>Line/Load Regulation</b>	±5%
<b>Leakage Current</b>	0.25mA max
<b>Inrush Current</b>	<60A @ 100Vac, 25°C , cold start <100A @ 240Vac, 25°C , cold start

### Environmental Specifications :

<b>Operating Temperature</b>	0°C ~ +40°C
<b>Storage Temperature</b>	-20°C ~ +75°C , non-condensing
<b>Operating Humidity</b>	20% ~ 90% RH
<b>Storage Humidity</b>	20% ~ 95% RH, non-condensing

### Safety and Compliance :

UL/cUL60950-1, SAAAS/NZS 60950-1:2003, CE approval
FCC Class B, EN55022 Class B, C-Tick approval
ESD : E61000-4-2, Level 4
Radiated Immunity: EN61000-4-3, Level 2
EFT: EN61000-4-4, Level 2
Surge: EN61000-4-5, Level 3
Conducted Immunity: EN61000-4-6, Level 2
Magnetic Immunity: EN61000-4-6, Level 1
Voltage Dips: EN61000-4-11

Note:

- Measurements shall be made with an oscilloscope with 20MHz bandwidth. Output shall be bypassed at a connector with a 0.1uF ceramic capacitor and a 10uF low ESR electrolytic capacitor.
- Efficiency: Measure efficiency at 100%, 75%, 50%, 25% load after burn-in 30 minutes at 100% load, then compute the arithmetic average of these 4 values.
- Interchangeable AC clip: US, Brazil, CN, EU, Korea, UK, India, Argentina, Australia, IEC320 C8

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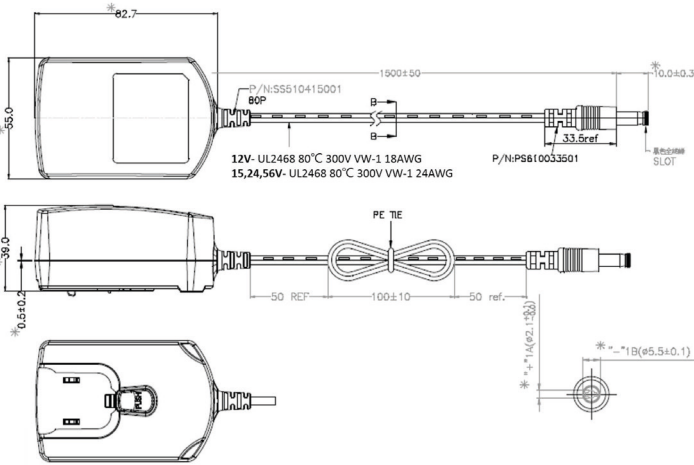
Source: www.phihong.com.tw, issue: 12.06.2017. Subject to change without notice.

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declares that the following product

Product name: **PROFILER SCD**  
 Article no.: **00-76-900-4830**  
 Product type: **Digital readout unit**  
**for 2D and 3D measurements**

is in conformity with the following European Union harmonisation legislation, when operated in accordance with the operating manual:

**EMC directive 2014/30/EU**  
**Low voltage directive 2014/35/EU**  
**RoHS directive 2011/65/EU**

The following harmonised standards have been applied:

**DIN EN 61010-1:2020-03**  
**DIN EN 61326-1:2013-07**

In case of a modification of the product, which was not approved by the manufacturer, this Declaration of Conformity loses its validity.

This Declaration of Conformity is issued under the sole responsibility of the manufacturer.

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Approved by Uwe Schnitter | 26.04.2022

Issued by Christopher Koch | 26.04.2022

Proved to be QMS conform by Uwe Schnitter | 26.04.2022

QM document AB502\_QF07\_07 | Rev. A | Ver. 00 | Page 1 of 1

## 16 Index

### A

accessories 112  
 acoustic signal 153  
 area of application 113  
 axis assignment 166

### B

baud rate 130, 167

### C

cables 112, 189  
 cleaning 181  
 COM  
   *button* 168  
   *menu* 167  
   *port* 171  
 configuration 152  
 conversion values 168  
 counting direction 164  
 Customer Service Center 188

### D

decimal places 135, 160  
 decl. of conformity 114, 204  
 display  
   *brightness* 154  
   *rotation* 160  
 Display (menu) 160  
 disposal 191  
 driver 126, 128

### E

encoder cables 112  
 encoder interface 194  
 Encoder (menu) 163  
 encoder type 164, 192  
 entry bar 134  
 evaluation options 163

### F

factory setting 169  
 file format 171, 179  
 firmware version 170  
 foot switch 131

### G

GetPos 171  
   *configuration* 173  
   *file format* 171, 179  
   *installation* 171  
   *key layout* 178  
   *menu functions* 176  
   *position tables* 179  
   *user interface* 175

### I

identification plate 113, 188  
 Information (menu) 170  
 installation requirements 116  
 instruction set 180  
 intended use 116

### L

language 154, 177

### M

machine zero 154  
 maintenance 182  
 measurement function  
   *enable/disable* 164  
 measurements 137  
 measuring system  
   *3rd* 122  
   *appropriate* 116, 192  
 measuring unit 135, 160

**O**

operating properties 153  
 operation 133  
 Origin (menu) 154

**P**

password query 152, 154  
 PC 126, 171  
 period 165  
 ports 114, 195  
 position correction  
   *configuration* 141  
   *delete* 150  
   *enable/disable* 150, 164  
   *functioning* 141  
   *modify* 148  
   *PosCorr (menu)* 161  
   *pre-configured* 141  
 position display 134  
 position value 154  
   *send* 131, 138, 176, 178  
   *zero* 131, 138, 154  
 power consumption 193  
 power supply 202  
 product description 113  
 putting into service 120

**R**

reference mark 155  
 reference signal evaluation 137  
   *enable/disable* 164  
 reference switch 155  
 replacement components 189  
 reset 169  
 RS-232 connection 129

**S**

safety 116  
 scope of delivery 111  
 security functions 194

service 188  
 Service (menu) 168  
 Setup (menu) 152  
 shield 201  
 signal amplitude 168  
 signs  
   *input field* 134  
   *position display* 164  
 standby mode 151  
 status display  
   *general* 136  
   *position correction* 136  
   *reference signal* 137  
 supply voltage 193  
 support 188  
 System (menu) 153  
 system requirements 193

**T**

taking out of service 190  
 technical data 192  
 terminal 130, 138  
 troubleshooting 183

**U**

USB connection 126  
 user interface 133  
 user training 188

**V**

version 170

**W**

wall installation 112, 160

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An den Fichten 35  
35579 Wetzlar (Germany)

Tel.: +49 6441 67106-0  
Fax: +49 6441 67106-28

[info@marzhauser-st.com](mailto:info@marzhauser-st.com)  
[www.marzhauser-st.com](http://www.marzhauser-st.com)

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© 2022 | *All rights reserved. Any complete or partial copy requires written authorization from Märzhäuser Sensotech GmbH.*
**Märzhäuser Sensotech GmbH**