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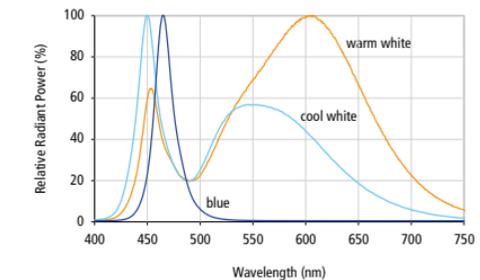
LED Illumination for Microscopy

Technical Data


Order no.	XX-59-11X-6800 cool white: 114 lm, 5,000–10,000 K
XX-59-11X-6810	warm white: 87.4 lm, 2,600–3,700 K
XX-59-11X-6820,	blue: 39.8 lm, 470–480 nm
16-59-110-6800	
XX: customer specific adapter, min. lumen at 350 mA	
Brightness	0.5–100 %, infinitely adjustable
Lumen maintenance	> 70 % after 20,000 hours, typ. (with sufficient cooling)
Angle of radiation	approx. 5°
Pulse interval	n*40 µs (TANGO + AUX I/O or DSU-1 + ext. trigger)
Operating voltage	10.8 ... 13.2 V
Protection type	IP43
Material	aluminium, black anodised
Dimensions	Ø 70 × 35 mm (without adapter)
Weight	approx. 0.5 kg

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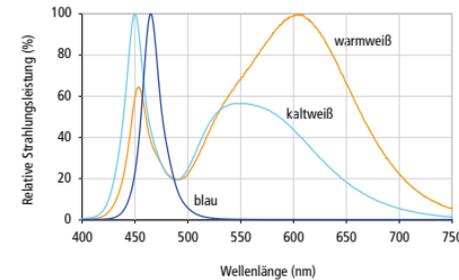Troubleshooting

The LED 100 does not work if e.g.

- *Digital OUT 0* is enabled in the *SwitchBoard* software.
- the DSU-1 is set to trigger mode.
- the connected DSU-1/TANGO is not powered.
- the dial control of the DSU-1 is set to *min*.

Service

You can reach our Customer Service Center Mondays to Fridays, from 8 am to 4 pm CET. Please keep handy the information stated on the type plate.

Tel.: +49 6441 9116-36

E-mail: service@marzhauser.com

Disposal

Referring to EU directive 2012/19/EU (WEEE) the LED 100 is classified as electronic equipment and may not be disposed of with household waste. Märzhäuser Wetzlar will properly dispose of the LED 100. Also note the regulations outside the EU.

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Declaration of Conformity

The LED 100 was designed and produced in compliance with applicable EU directives and harmonized standards. The Declaration of Conformity is available on request at:

service@marzhauser.com

Cleaning

WARNING

Danger to life from spreading hazardous substances.

In case of possible contamination with corrosive, infectious, toxic, radioactive or other harmful substances disinfect/decontaminate the LED 100 professionally.

CAUTION

Risk of burns when touching the LED 100 after long operation time.

Let the LED 100 cool down.

NOTICE

Risk of equipment damage from improper cleaning.

Ensure that no liquids enter the LED 100 or its plug.

1 Take the LED 100 out of operation.

2 Clean the LED 100 with a lint-free cleaning cloth. With greater contamination, slightly dampen the cleaning cloth with water or a mild solvent-free disinfectant/cleaning agent.

3 Allow the LED 100 to dry.

Taking Out of Service

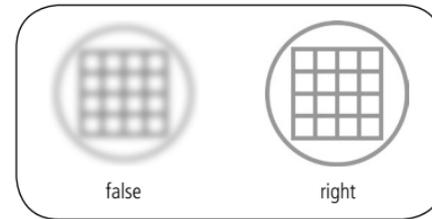
1 Switch off the DSU-1/TANGO and the microscope.

2 Disconnect the LED 100 from the DSU-1/TANGO.


3 Unmount the LED 100 from the microscope.

4 Clean the LED 100.


5 Switch on the DSU-1/TANGO and evaluate the image.

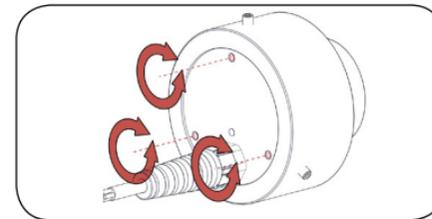6 Adjust the LED 100 using the 3 screws on the back until the image matches with the figure „right“. Make sure that the back of the LED 100 sits in the locking points (every 120°) of the housing.

7 Tighten the 3 screws.

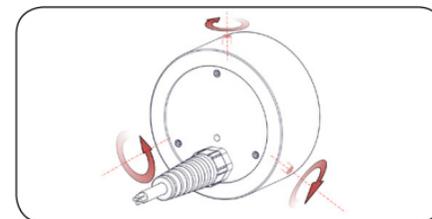8 Switch off the DSU-1/TANGO.

9 Mount the LED 100 to the microscope.

Putting into Service

1 Switch off the DSU-1/TANGO and the microscope.

2 Connect the LED 100 with the *LED 100* port of the DSU-1 or the *AUX I/O* port of the TANGO.

3 Put the DSU-1/TANGO into operation as described in the respective operating manual.

Adjustment (optional)

In exceptional cases, such as a weak illumination of the microscope, the LED 100 has to be adjusted.

CAUTION

Risk of eye damage from very bright light.

Do not look into the LED 100 during operation.

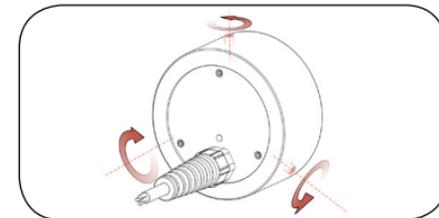
Risk of burns when touching the LED 100 after long operation time.

Let the LED 100 cool down.

1 Switch off the DSU-1/TANGO and the microscope.

2 Unmount the LED 100 from the microscope.

3 Hold the housing and the back of the LED 100 tightly with your hand. Loosen the 3 screws.

4 Direct the LED 100 towards a plain surface (e.g. white wall) in a distance of approx. 5 meters.

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General Safety Instructions

The LED illumination LED 100 is intended for the illumination of a microscope and must be operated with a Märzhäuser controller (DSU-1 or TANGO) only.

- Read and understand this operating manual as well as the operating manual of the DSU-1/TANGO and the microscope.
- The LED 100 is in accordance with DIN EN 62471:2009 „Photobiological Safety of Lamps“ classified in Risk Group 2. Do not look directly into the light of the LED 100 in order to avoid eye damage. Using the LED 100 on a microscope is not dangerous as long as the aversion response to bright light is given. Decrease the intensity of the LED 100 as much as necessary, so that you are not blinded during microscoping.
- In trigger mode, pulsing/flashing light can cause epileptic reactions.
- Disconnect the DSU-1/TANGO from the mains before connecting/disconnecting the LED 100.
- Overheating of the LED 100 can have a negative effect on the service life. Therefore, do not encase the LED 100 and do not cover the housing of the LED 100.
- Ensure that no chips, liquids, objects or substances enter the interior of the LED 100. Protect the LED 100 from dust and dirt as much as possible.
- The LED 100 is not protected against corrosive, infectious, toxic, radioactive or other substances, which are harmful to health. When handling such substances, ensure that all legal requirements are fulfilled, especially the national regulations on accident prevention.
- Immediately replace the LED 100 if there is damage on the housing or cable.
- Open the housing of the LED 100 for adjusting reasons only.

- Only have repair work carried out by our Customer Service Center (see chapter *Service*).
- When lending or selling the LED 100, make sure to provide all the documents supplied with it.

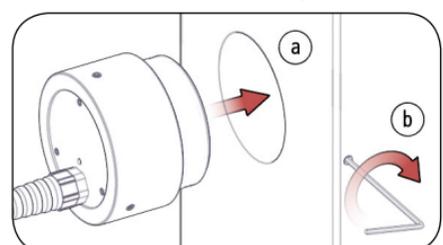
Delivery Contents

LED 100, hex key, operating manual

Assembly

 CAUTION
 Risk of eye damage from very bright light. Do not look into the LED 100 during operation.
 Risk of burns when touching the LED 100 after long operation time. Let the LED 100 cool down.

- 1 Switch off the microscope.
- 2 If necessary, remove the lamp housing of the microscope as described in the operating manual.
- 3 Mount the LED 100 to the microscope.

Note: Depending on the adapter the assembly of the LED 100 can be different from the figure.

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Märzhäuser Wetzlar GmbH & Co. KG