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Operating Manual DSU-1

Controller for LED 100

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12 Watt Interchangeable Plug Series

Features

- Double Insulated
- Class B EMI
- Level VI Efficiency Compliant
- Limited Power Source (LPS)

Applications

- Wireless Communications
- Network Equipment
- Peripherals
- Gaming

Safety Approvals

- cUL/UL 60950-1
- cUL/UL 62368-1
- AS/NZ 60950-1
- IEC60950-1
- IEC62368-1
- CE

Mechanical Characteristics

- Length: 71.6mm (2.82in)
- Width: 50mm (1.97in)
- Height: 33.25mm (1.22in)
- Weight: 112g (3.95oz)

Output Specifications

Model	DC Output Voltage	Load		Ripple ⁽¹⁾ P-P (max.)
		Min.	Max.	
PSAC12R-050-R	5V	0A	2.4A	100mV
PSAC12R-060-R	5.9V	0A	2.0A	100mV
PSAC12R-075-R	7.5V	0A	1.6A	100mV
PSAC12R-090-R	9V	0A	1.33A	120mV
PSAC12R-120-R	12V	0A	1.0A	120mV
PSAC12R-150-R	15V	0A	0.8A	150mV
PSAC12R-240-R	24V	0A	0.5A	240mV
PSAC12R-480-R	48V	0A	0.25A	480mV

Notes:

1. Ripple measured by using a 12-inch twisted pair terminated with 10uF capacitor and 0.1uF ceramic in parallel with oscilloscope set to 20Mhz measured after a warm-up of 10minutes.

Phihong is not responsible for any error and reserves the right to make changes without notice. Please visit our website at www.phihong.com for the most up-to-date specifications and contact information. Revision 1/16/2020

PSAC12R Characteristics ¹		WWW.PHIHONG.COM
INPUT		Immunity
Input Voltage Rating	100 to 240VAC	EN61000-4-2
		EN61000-4-3
		EN61000-4-4
AC Input Voltage Range	90 to 264VAC	EN61000-4-5 Level 3
		EN61000-4-6
		EN61000-4-11
AC Input Current	0.5A (RMS), 120VAC at maximum load	Isolation (HI-POT test)
	0.25A (RMS), 230VAC at maximum load	Input to Output: 3000VAC for 1 minute, 10mA
AC Input Frequency	47 to 63Hz	Insulation Resistance
		Input to Output: 500VDC 500M Ohm minimum
In-rush Current	<30A for 120VAC at maximum load	Over Current Protection
	<60A for 230VAC at maximum load	3A max. Short Circuit Auto-restart
	(cold start at ambient 25°C)	Short Circuit Protection
		Output can be shorted without damage
Leakage Current	0.25mA maximum	Over Voltage Protection
		<7.5V
OUTPUT		Dielectric Withstand (Hi-pot) Test
Output Power	12W	Primary to Secondary: 3000V AC for 1 min., 10mA
Efficiency²		DC Cord
US DoE Level VI		1500mm (18~24AWG)
EU CoC Ver5 Tier1		DC Output Connector (Tuning Fork Type)
No Load Power Consumption	≤75mW max @ 115V	5.5mm x 2.1mm x 10mm Center Positive
Hold up time	10mS minimum at maximum load, 120VAC	Interchangeable AC Clips
		(Sold Separately)
ENVIRONMENTAL		RPA – US
Temperature		RPB – Brazil
Operation	0°C to +40°C	RPC – China
Non-operation	-40°C to +85°C	RPE – Europe
Humidity	50 to 90%	RPH – Korea
		RPI – India
		RPK – UK
EMI		RPN – Argentina
Complies with FCC Class B		RPS – Australia
Complies with EN55032 Class B		RPX – IEC320 C8

Notes:

1. The characteristics defined are at ambient temperature of 25°C unless otherwise specified
2. Efficiency is measured after 30 minutes burn-in

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hereby declares that the following product
 Product name: **DSU-1**
 Product type: **Controller**
 Indented use: **Dimmer-Shutter Unit**
for LED 100
 Item no.: **00-76-110-6841**

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is in conformity with the following relevant
 European Union harmonisation legislation:
 until 19.04.2016
EMC Directive 2004/108/EC
Low Voltage Directive 2006/95/EC
RoHS Directive 2011/65/EU
 from 20.04.2016
EMC Directive 2014/30/EU
Low Voltage Directive 2014/35/EU
RoHS Directive 2011/65/EU

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This declaration of conformity is issued under the
 sole responsibility of the manufacturer.
 In case of a modification of the product, which was
 not approved by the manufacturer, this declaration
 loses its validity.

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Approval:	QMS-Conform:	Issued by:	Time:
16:20 Date	16:20 Date	Schnitter Uwe Date / Sign:	16:20
 Signature	 Signature	16:20 / sch	

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Wartung 23
WEEE 29

Table of Contents

1	Basics of the Operating Manual	39
1.1	Symbols and Abbreviations	40
2	Delivery Contents	41
3	Product Description	42
3.1	Identification	42
3.2	Area of Application	42
3.3	Compliance with Directives and Standards	43
3.4	Structure and Ports	44
3.5	Trigger Function (Shutter)	46
4	Safety	47
4.1	Intended Use	47
4.2	Installation Requirements	47
4.3	General Safety Instructions	48
4.4	Warning Note Categories	50
4.5	Used Hazard Symbols	50
5	Commissioning	51
6	Operation	53
7	Cleaning	54
8	Maintenance	55
9	Troubleshooting	55
10	Service	58
10.1	Replacement Components	59
11	Decommissioning	60
12	Disposal	61
13	Technical Data	62
13.1	DSU-1	62
13.2	External 12 V Power Supply (PSAC12R-120-R)	64
14	EC Declaration of Conformity	66
15	Index	67

1 Basics of the Operating Manual


This operating manual gives you the information which is needed for trouble-free and safe use of the DSU-1 (Dimmer-Shutter Unit). You must have fully read and understood this operating manual and especially the safety instructions before you commission, operate, clean, decommission or dispose of the DSU-1.

Märzhäuser Sensotech accepts no liability for damage and operating faults which result from ignoring this operating manual.

1.1 Symbols and Abbreviations

Symbol	Meaning
	additional information and useful notes
	start of an instruction
1, 2, 3 ...	steps within an instruction
»	cross reference to another part of this operating manual, another document or a website
Abbreviation	Meaning
approx.	approximately, roughly
e.g.	for example
fig.	figure
LED	light-emitting diode
max.	maximum
min.	minimum
resp.	respectively
typ.	typical

2 Delivery Contents

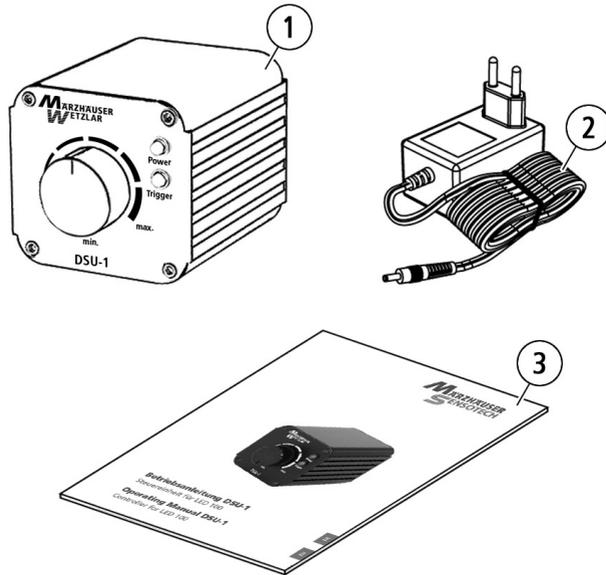No.	Designation	Quantity
1	DSU-1 (Dimmer-Shutter Unit)	1
2	External 12 V power supply incl. 4 adapters (EU/GB/USA/AUS, no fig.)	1
3	Operating manual	1

Fig. 2: © Michael E. Kimbler/Markus Walter, source: 3D Content Central, Dassault Systèmes

3 Product Description

3.1 Identification

You can find all required information for clear identification of the DSU-1 on the identification plate. The identification plate is attached to the underside of the DSU-1.

3.2 Area of Application

The DSU-1 is a control unit for the LED illumination LED 100. With the DSU-1, the LED 100 can be operated independently and without a TANGO controller.

By using the dial control the light intensity is adjustable from 0.5...100 %. A digital trigger input also enables quick switching of the LED 100 and the realization of a trigger function (shutter).

The DSU-1 is powered by an external 12 V power supply which also powers the LED 100.

Here you can find further information on the LED 100:

» www.marzhauser.com/led100

3.3 Compliance with Directives and Standards

 Märzhäuser Sensotech hereby declares that the DSU-1 was designed and produced in compliance with the EC/EU directives and harmonized standards mentioned in the EC Declaration of Conformity.

The EC Declaration of Conformity can be found here:

» *Chapter 14 EC Declaration of Conformity, page 66*

3.4 Structure and Ports

Front panel

Rear panel

No.	Designation	Function
1	Dial control	adjust light intensity from 0.5...100 %
2	Power	LED is on when DSU-1 is switched on
3	Trigger	LED is on when trigger function (shutter) is switched on
4	Power ON/OFF	on-off switch ¹
5	Trigger ON/OFF	on-off switch for trigger function (shutter)
6	12V / 0,5A max.	port for power supply
7	Trigger IN	port for external trigger signal generator
8	LED 100	port for LED 100

¹ **Please note:** Switching off the DSU-1 causes a disconnection of the LED 100 from the mains. To completely disconnect the DSU-1 from the mains, remove the power supply from the mains.

3.5 Trigger Function (Shutter)

The trigger function (shutter) enables a rapid on-off switching of the LED 100 by means of a digital control signal. Due to the short on-off times and the very low afterglow time of the light source, a precise shutter function can be realized.

The digital control signal can be connected to the *Trigger IN* signal input port. A high signal at the signal input port turns off the LED 100. If a low signal is transmitted or if no external trigger signal generator is connected, the LED 100 shines according to the selected light intensity. During the trigger process, the light intensity of the LED 100 can still be controlled by means of the dial control.

The trigger function can be enabled/disabled by means of the *Trigger ON/OFF* switch. When the trigger function is enabled, the *Trigger LED* is on.

Technical specifications of the trigger function can be found here:

» *Chapter 13 Technical Data, page 62*

4 Safety

4.1 Intended Use

The DSU-1 (Dimmer-Shutter Unit) is intended and designed as an independent

- power supply and
- control unit

for the LED illumination LED 100.

4.2 Installation Requirements

Category	Condition
Installation location	closed rooms, no risk of explosion
Installation height	max. 2,000 m
Environmental temperature	0...+40 °C
Storage temperature	-20...+70 °C
Cooling	normal convection
Humidity	max. 85 %, non-condensing
Protection type	IP40

4.3 General Safety Instructions

- Only use the DSU-1 for the purposes which are described in this operating manual.
- Read and understand the operating manuals of the LED 100 and peripherals, which you would like to connect to the DSU-1.
- Only connect light sources type LED 100 from Märzhäuser Wetzlar to the *LED 100* port.
- Only use shielded cables for signal transfer.
- Only use the supplied external 12 V power supply to power the DSU-1.
- Ensure that the mains voltage matches the label on the power supply before connecting the DSU-1 to the mains.
- Disconnect the DSU-1 from the mains before connecting/ disconnecting the LED 100 or an external trigger signal generator.
- Set the DSU-1 in such a way that you can immediately disconnect the power supply from the mains, without hindrance, in case of emergency.
- Ensure that no chips, liquids, objects or substances enter the interior of the DSU-1 or the power supply. Protect the DSU-1 from dust and dirt as much as possible.
- DSU-1 is not protected against corrosive, infectious, radioactive or other substances, which are harmful to health. When handling such substances, ensure that all legal requirements are fulfilled, especially the national regulations on accident prevention.
- Immediately replace the power supply if there is damage on the housing or cable of the power supply.
- Disconnect the DSU-1 from the mains if there are malfunctions, defects or safety flaws. Secure the DSU-1 against restarting.

- Never open the housing of the DSU-1 or the power supply. In case of violation, all warranty and liability claims will become void.
- Only have repair work carried out by our Customer Service Center.
 - » *Chapter 10 Service, page 58*
- Ensure that the DSU-1, in conjunction with your application, meets the applicable safety regulations and legal requirements.
- If you lend out or sell the DSU-1, provide all the documents supplied with it.

4.4 Warning Note Categories

Warning notes are divided into the following categories:

WARNING

Hazardous situation that can cause serious injury or death if not avoided

NOTICE

Warning of possible damage or malfunction

4.5 Used Hazard Symbols

The following hazard symbols are used in warning notes:

Symbol	Meaning
	warning of hazard from electric current
	warning of hazard from substances which are harmful to health

5 Commissioning

WARNING

Possible electric shock and equipment damage from connecting non-specified external trigger signal generators!

The DSU-1 is insulated according to the power supply. Ensure that no dangerous voltages or currents are induced in the DSU-1 by an external trigger signal generator.

Possible electric shock and equipment damage from connecting foreign or damaged power supplies!

Only use the supplied external 12 V power supply. Immediately replace the power supply if there is damage on the housing or cable of the power supply. Only use replacement components from Märzhäuser Sensotech.

Possible electric shock and equipment damage from connecting to non-specified mains!

Ensure that your mains voltage matches the label on the power supply before connecting the DSU-1 to the mains.

NOTICE

Possible equipment damage from commissioning/decommissioning with existing power connection!


Disconnect the DSU-1 from the mains before connecting/disconnecting the LED 100 or an external trigger signal generator.

Possible incorrect data transmission from using an unshielded cable!

Only use a shielded cable for connecting an external trigger signal generator.

► Commissioning the DSU-1

- 1 Ensure that the DSU-1 has a secure footing.
- 2 Ensure that the DSU-1 is switched off (*Power ON/OFF* switch in *OFF* position).
- 3 Set the dial control to position *min.*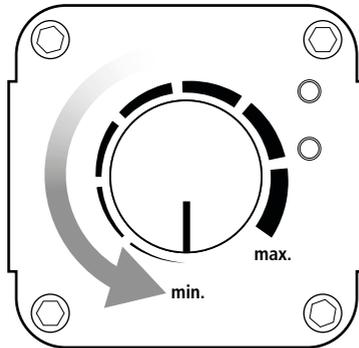

- 4 Connect the cable of the LED 100 to the *LED 100* port.
- 5 If needed, connect the cable of the external trigger signal generator to the *Trigger IN* port.
- 6 Connect the power supply cable to the *12V / 0,5A max.* port.
- 7 Connect the power supply to the mains.

The *Power* LED lights up and signals that the DSU-1 is ready for operation. Commissioning is now complete.

6 Operation

► Switching on/off the DSU-1

To switch on the DSU-1, set the *Power ON/OFF* switch to the *ON* position. The *Power* LED lights up.

To switch off the DSU-1, set the *Power ON/OFF* switch to the *OFF* position. The *Power* LED turns off.

► Enabling/disabling the trigger function (shutter)


To enable the trigger function (shutter), set the *Trigger ON/OFF* switch to the *ON* position. The *Trigger* LED lights up.

To disable the trigger function (shutter), set the *Trigger ON/OFF* switch to the *OFF* position. The *Trigger* LED turns off.

► Controlling the light intensity

In order to increase the light intensity of the LED 100, turn the dial control clockwise.

In order to decrease the light intensity of the LED 100, turn the dial control counterclockwise.

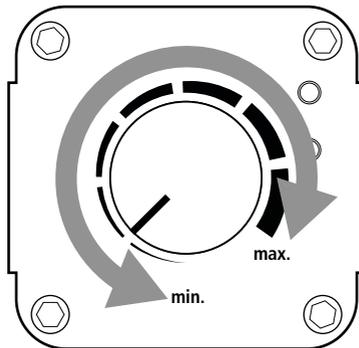7 Cleaning

WARNING

Danger to life and risk of equipment damage from improper cleaning!

Disconnect the power supply from the mains before cleaning the DSU-1 or the power supply. Ensure that no liquids enter the interior of the DSU-1 or the power supply. Protect ports and plugs from liquids. Never touch the power supply or the socket with wet hands.

Danger to life and risk of equipment damage from spreading hazardous substances!

In case of possible contamination with corrosive, infectious, toxic, radioactive or other substances which are harmful to health, have the DSU-1 professionally disinfected/decontaminated.

► **Cleaning the DSU-1**

- 1 Switch off the DSU-1.
- 2 Disconnect the power supply from the mains.
- 3 Clean the DSU-1 and the power supply with a lint-free cleaning cloth. With greater contamination, slightly dampen the cleaning cloth with water or a mild solvent-free disinfectant/cleaning agent.
- 4 Allow the DSU-1 and the power supply to dry.
- 5 Connect the power supply to the mains.
- 6 Switch on the DSU-1.

The DSU-1 can now be used again.

8 Maintenance

The DSU-1 is maintenance-free.

9 Troubleshooting

WARNING

Possible electric shock and equipment damage from connecting a foreign or damaged power supply!

Only use the supplied external 12 V power supply. Immediately replace the power supply if there is damage on the housing or cable of the power supply. Only use replacement components from Märzhäuser Sensotech.

Possible electric shock and equipment damage from not permitted repair attempts!

Never open the housing of the DSU-1 or the power supply. Only have repair work carried out by our Customer Service Center. In case of violation, all warranty and liability claims will become void.

The following overview can help you in removing faults and malfunctions of the DSU-1.

If a fault or malfunction cannot be remedied using the overview, contact our Customer Service Center.

» *Chapter 10 Service, page 58*

Fault	Possible causes
DSU-1 can not be switched on	<p data-bbox="594 229 706 256">no power</p> <hr/> <p data-bbox="594 432 1006 459">power supply or DSU-1 out of order</p>
Connected LED 100 does not work although <i>Power</i> LED is on	<p data-bbox="594 533 882 560">light intensity set to <i>min.</i></p> <hr/> <p data-bbox="594 635 837 662">LED 100 out of order</p> <hr/> <p data-bbox="594 735 1005 826">trigger function (shutter) enabled, signal generator sends a high signal continuously</p> <hr/> <p data-bbox="594 874 882 901">incompatible light source</p>
Limited or no trigger function (shutter)	<p data-bbox="594 1007 978 1034">trigger function (shutter) disabled</p> <hr/> <p data-bbox="594 1107 740 1134">faulty wiring</p>

Remedy

- 1 Make sure that the voltage of your mains corresponds to the label on the power supply.
 - 2 Make sure that the power supply is properly connected to the DSU-1 (12V / 0.5 A max. port) and the mains.
 - 3 Make sure that the needed mains voltage is available.
-

Contact our Customer Service Center.

» *Chapter 10 Service, page 58*

Check the light intensity setting.

» *Chapter 6 Operation, page 53*

Contact our Customer Service Center.

» *Chapter 10 Service, page 58*

Check the sended trigger signal or disable the trigger function (shutter).

» *Chapter 6 Operation, page 53*

» *Chapter 13.1 DSU-1, page 62*

Make sure that the light source corresponds to the specifications of the DSU-1 and is approved by Märzhäuser Sensotech.

» *Chapter 4.1 Intended Use, page 47*

Aktivate the trigger function (shutter).

» *Chapter 6 Operation, page 53*

- 1 Make sure that the cable is shielded.
 - 2 Make sure that the external trigger signal generator is working properly and connected to the *Trigger IN* port.
 - » *Chapter 5 Commissioning, page 51*
-

10 Service

The staff at our Customer Service Center are happily at your disposal for support queries, complaints and suggestions about our products and services.

Overview of the services from our Customer Service Center:

- Advice and support
- Support for customer-specific application questions
- Support for commissioning equipment
- Inspection and repair of equipment
- Error analysis (also online via web conference)
- Delivery of replacements or replacement components
- User training (also online via web conference)

You can reach our Customer Service Center Mondays to Fridays from 08:00 am to 04:00 pm CET.

Tel.: **+49 6441 9116-36**
 Fax: **+49 6441 9116-40**
 E-mail: **service@marzhauser.com**

So that we can help you quickly when you need service, please keep handy the following information about your DSU-1:

Description	Information on the identification plate
Type	
S/N	
Art.-No.	

10.1 Replacement Components

Should you need to replace a component of the DSU-1, please contact our Customer Service Center.

» *Chapter 10 Service, page 58*

Replacement components	Order no.
External 12 V power supply	00-76-110-6842
Operating manual	00-76-110-6841-TD

11 Decommissioning

► Decommissioning the DSU-1

- 1 Switch off the DSU-1.

NOTICE

Risk of equipment damage from commissioning/decommissioning while the DSU-1 is connected to the mains!

Disconnect the DSU-1 from the power supply before connecting/disconnecting the LED 100 or an external trigger signal generator.

-
- 2 Disconnect the power supply of the DSU-1 from the mains.
 - 3 Disconnect all cables from the DSU-1.
 - 4 Clean the DSU-1.
 - » *Chapter 7 Cleaning, page 54*

The DSU-1 can now be disposed of.

12 Disposal

The adjacent symbol means that the DSU-1 is classified as electrical or electronic equipment according to EU directive 2012/19/EU (WEEE). The DSU-1, therefore, may not be disposed of together with household waste.

Since the DSU-1 is a commercially used device, disposal at municipal collection points for electrical and electronic equipment is prohibited.

Send the DSU-1 to Märzhäuser Sensotech for cost-free and proper disposal.

► Disposing of the DSU-1

- 1 Decommission the DSU-1.
» *Chapter 11 Decommissioning, page 60*
- 2 Send the DSU-1 for disposal to:
Märzhäuser Sensotech GmbH
An den Fichten 35
35579 Wetzlar (Germany)

All equipment purchased from Märzhäuser Sensotech can always be sent back for free disposal. If you dispose of a device outside the EU, note and follow the respective legal provisions.

13 Technical Data

13.1 DSU-1

Features

Control of light intensity	0.5...100 %
Dial control	industrial potentiometer with positive logarithmic characteristic
Operating voltage	11.4...12.6 V
Current consumption	$U = 12 \text{ V}, I_E = 0.4 \text{ A}$
Shielding	shield on housing
Protection type	IP40
Housing	aluminum
Dimensions	108 × 66 × 71 mm (L × W × H)
Weight	approx. 400 g (without accessories)

Trigger function (Shutter)

Port	Trigger IN
Input signal	TTL
Polarity	Active low

Trigger function (Shutter)

Turn on threshold	LED off: $U \geq 2,4 \text{ V}$ LED on: $U \leq 0,8 \text{ V}$ or open
-------------------	---

Internal resistance	$R_i \geq 10 \text{ k}\Omega$
---------------------	-------------------------------

Switch-on delay	approx. $35 \mu\text{s}$
-----------------	--------------------------

Switch-off delay	approx. $25 \mu\text{s}$
------------------	--------------------------

If needed, we provide measuring reports to support you using the trigger function. Don't hesitate to contact our Customer Service Center.

» *Chapter 10 Service, page 58*

13.2 External 12 V Power Supply (PSAC12R-120-R)

WWW.PHIHONG.COM

12 Watt Interchangeable Plug Series

Features

- Double Insulated
- Class B EMI
- Level VI Efficiency Compliant
- Limited Power Source (LPS)

Applications

- Wireless Communications
- Network Equipment
- Peripherals
- Gaming

Safety Approvals

- cUL/UL 60950-1
- cUL/UL 62368-1
- AS/NZ 60950-1
- IEC60950-1
- IEC62368-1
- CE

Mechanical Characteristics

- Length: 71.6mm (2.82in)
- Width: 50mm (1.97in)
- Height: 33.25mm (1.22in)
- Weight: 112g (3.95oz)

Output Specifications

Model	DC Output Voltage	Load		Ripple ⁽¹⁾ P-P (max.)
		Min.	Max.	
PSAC12R-050-R	5V	0A	2.4A	100mV
PSAC12R-060-R	5.9V	0A	2.0A	100mV
PSAC12R-075-R	7.5V	0A	1.6A	100mV
PSAC12R-090-R	9V	0A	1.33A	120mV
PSAC12R-120-R	12V	0A	1.0A	120mV
PSAC12R-150-R	15V	0A	0.8A	150mV
PSAC12R-240-R	24V	0A	0.5A	240mV
PSAC12R-480-R	48V	0A	0.25A	480mV

Notes:

1. Ripple measured by using a 12-inch twisted pair terminated with 10uF capacitor and 0.1uF ceramic in parallel with oscilloscope set to 20Mhz measured after a warm-up of 10minutes.

Phihong is not responsible for any error and reserves the right to make changes without notice. Please visit our website at www.phihong.com for the most up-to-date specifications and contact information. Revision 1/16/2020

PSAC12R Characteristics ¹		WWW.PHIHONG.COM
INPUT		Immunity
Input Voltage Rating	100 to 240VAC	EN61000-4-2 EN61000-4-3 EN61000-4-4 EN61000-4-5 Level 3 EN61000-4-6 EN61000-4-11
AC Input Voltage Range	90 to 264VAC	
AC Input Current	0.5A (RMS), 120VAC at maximum load 0.25A (RMS), 230VAC at maximum load	Isolation (HI-POT test) Input to Output: 3000VAC for 1 minute, 10mA
AC Input Frequency	47 to 63Hz	Insulation Resistance Input to Output: 500VDC 500M Ohm minimum
In-rush Current	<30A for 120VAC at maximum load <60A for 230VAC at maximum load (cold start at ambient 25°C)	Over Current Protection 3A max. Short Circuit Auto-restart
Leakage Current	0.25mA maximum	Short Circuit Protection Output can be shorted without damage
OUTPUT	Output Power 12W	Over Voltage Protection <7.5V
Efficiency²	US DoE Level VI EU CoC Ver5 Tier1	Dielectric Withstand (Hi-pot) Test Primary to Secondary: 3000V AC for 1 min., 10mA
No Load Power Consumption	≤75mW max @ 115V	DC Cord 1500mm (18~24AWG)
Hold up time	10mS minimum at maximum load, 120VAC	DC Output Connector (Tuning Fork Type) 5.5mm x 2.1mm x 10mm Center Positive
ENVIRONMENTAL		Interchangeable AC Clips (Sold Separately) RPA – US RPB – Brazil RPC – China RPE – Europe RPH – Korea RPI – India RPK – UK RPN – Argentina RPS – Australia RPX – IEC320 C8
Temperature	Operation 0°C to +40°C Non-operation -40°C to +85°C Humidity 50 to 90%	
EMI	Complies with FCC Class B Complies with EN55032 Class B	

Notes:

1. The characteristics defined are at ambient temperature of 25°C unless otherwise specified
2. Efficiency is measured after 30 minutes burn-in

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© 2020 | *All rights reserved. Any complete or partial copy requires written authorization from Märzhäuser Sensotech GmbH.*